


Cadenza CW-C18

150 x 2 mm

Application

## LC-MS/MS Application for Sphingolipids (Sphinganine-1-PO<sub>4</sub>)

スフィンゴ脂質(スフィンガニン-1-リン酸)のLC-MS/MS分析

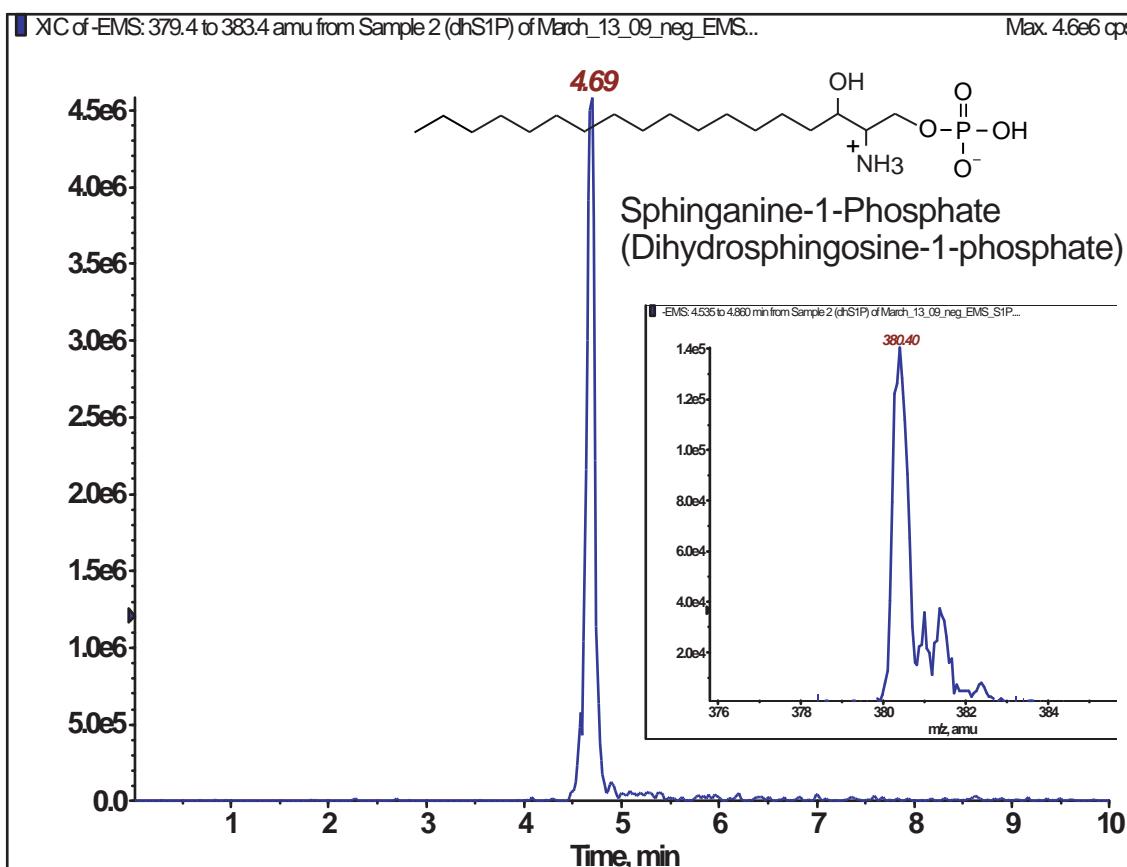

Cadenza CW-C18, 150 x 2 mm

A: [water / HCOOH (100:0.2 v/v)] containing 7mM ammonium formate  
 B: [methanol / HCOOH (100:0.2 v/v)] containing 5mM ammonium formate  
 70-90%B (0-1min), 90%B (1-1.5min), 90-100%B (1.5-2.5min),  
 100%B (2.5-6min), 100-40%B (6-6.5min), 40%B (6.5-10min)  
 0.2mL/min - 0.28mL/min (0-6min), 0.2mL/min (6-10min)  
 48deg.C, 5uL (50ug / mL)

Applied Biosystems QTrap, ESI Negative

Courtesy of Jennifer Ro and Joe Williams  
 Department of Evolution, Ecology, and Organismal Biology, Ohio State University - USA