



Unison UK-C8

100 x 3 mm

Application

LC-MS: low-calorie sweeteners

低カロリー甘味料のLC-MS分析

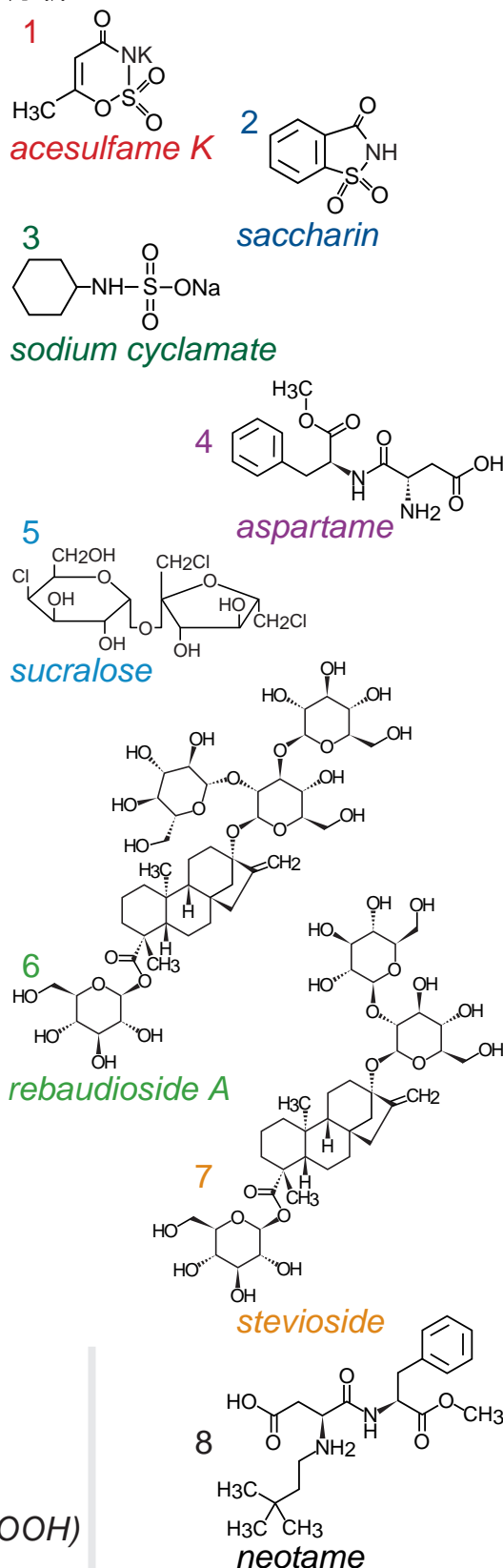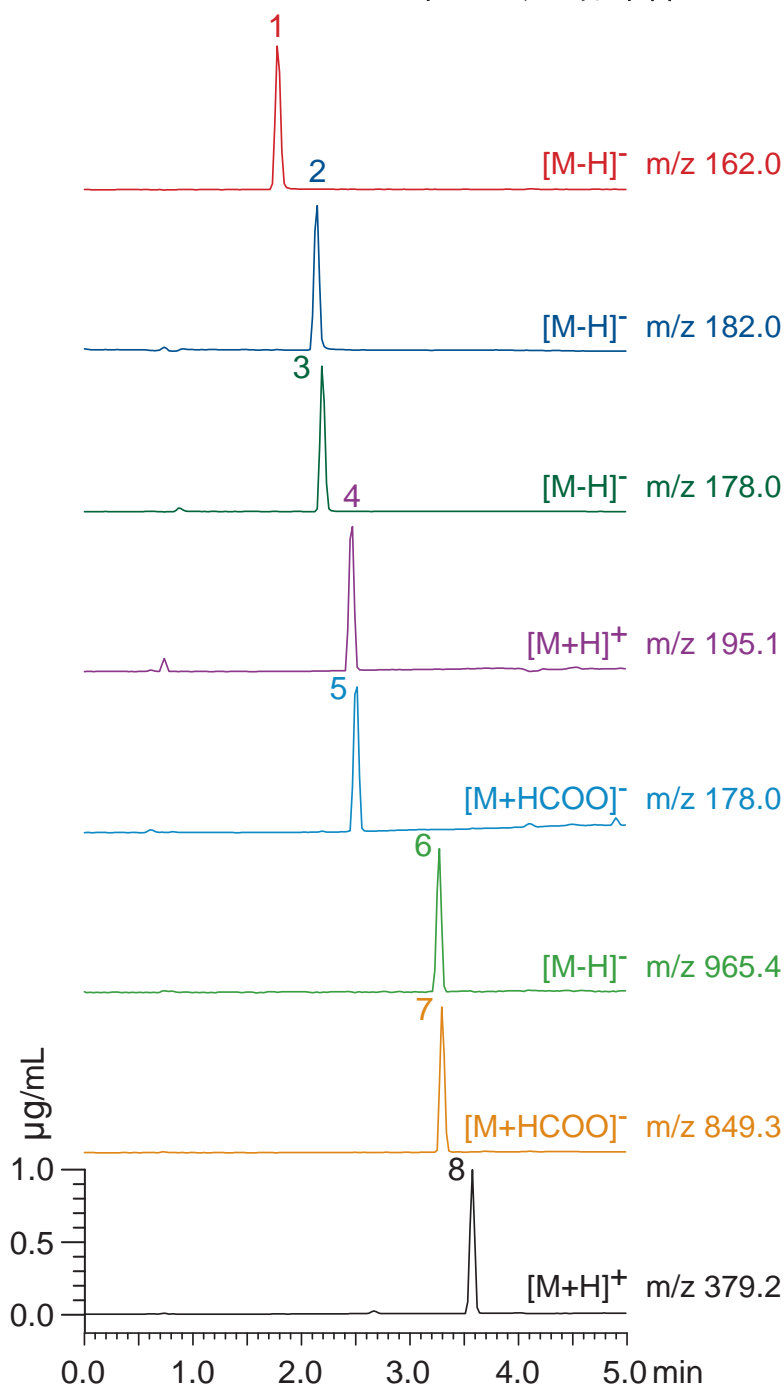

Unison UK-C8, 100 x 3 mm
 A: water / formic acid = 100 / 0.1
 B: acetonitrile / formic acid = 100 / 0.1
 10-80 %B (0-5 min), 100 %B (5-6 min)
 0.6 mL/min (10 MPa), 37°C, 2 µL(1µg/mL, 0.1% HCOOH)
 Single Quad. MS (ESI, positive, negative)