




Intrada Amino Acid
Scherzo SM-C18

50 x 3 mm

Application

LC-MS: Thyroid hormone


甲状腺ホルモン

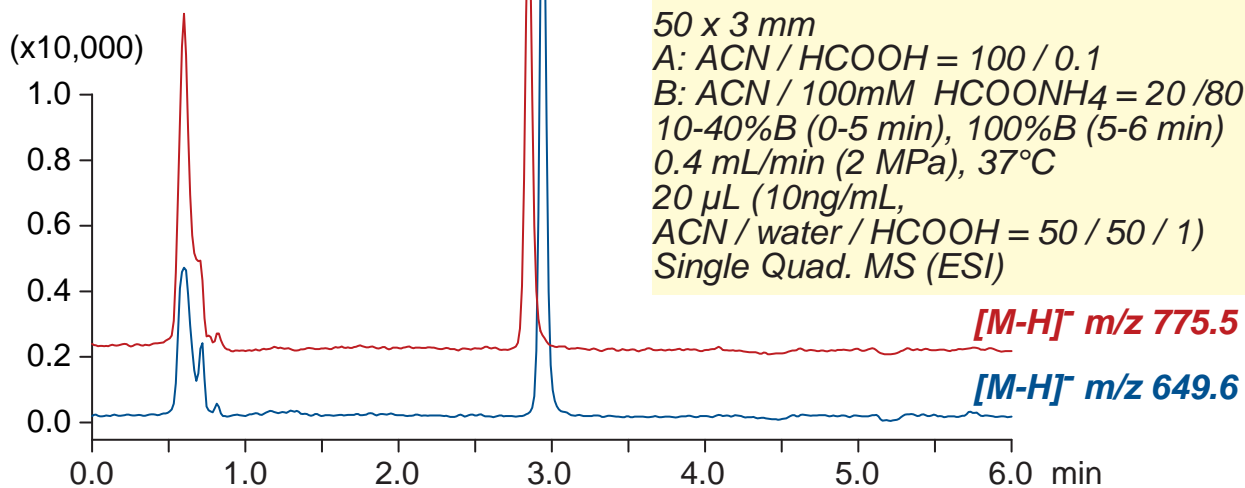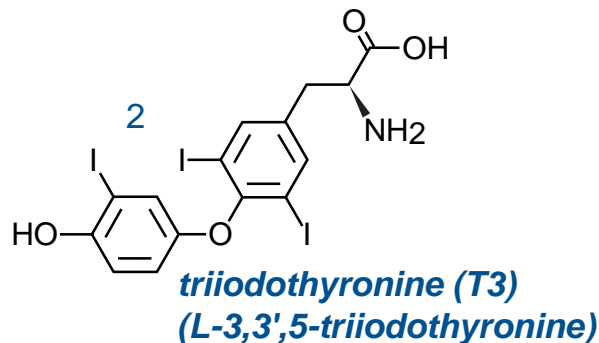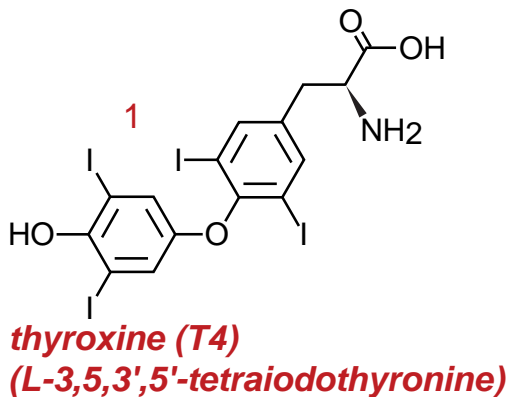

Intrada Amino Acid

50 x 3 mm
A: ACN / HCOOH = 100 / 0.1
B: ACN / 100mM HCOONH₄ = 20 / 80
10-40%B (0-5 min), 100%B (5-6 min)
0.4 mL/min (2 MPa), 37°C
20 μL (10ng/mL,
ACN / water / HCOOH = 50 / 50 / 1)
Single Quad. MS (ESI)

Scherzo SM-C18

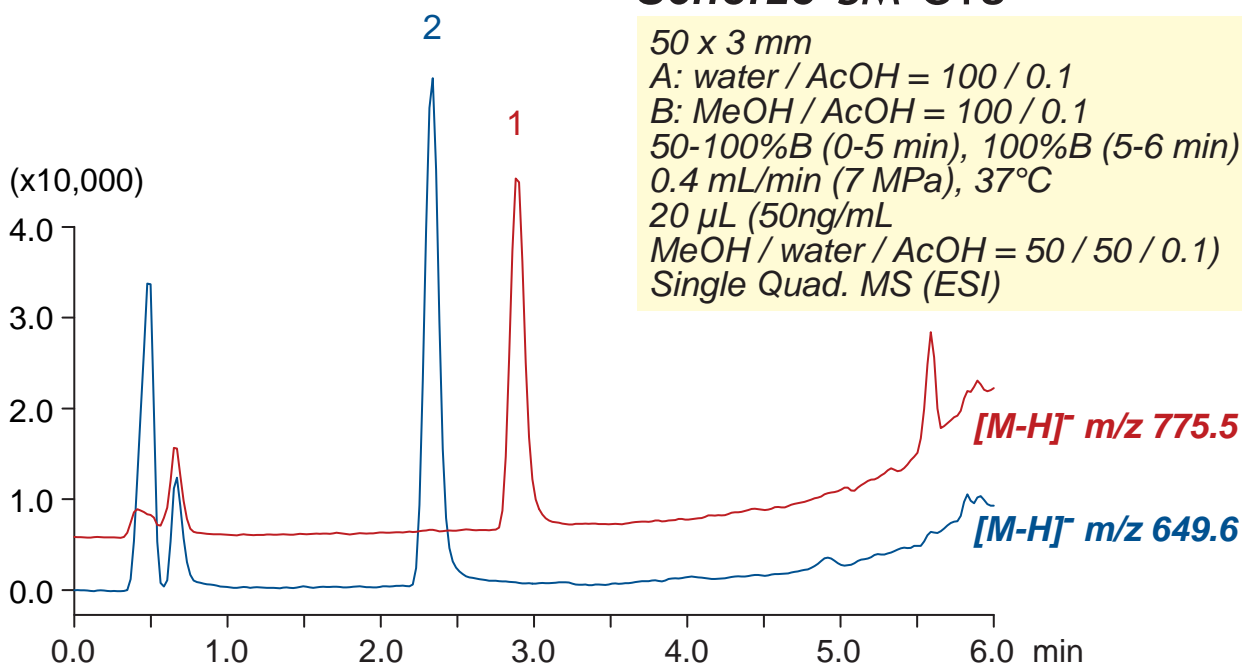

50 x 3 mm
A: water / AcOH = 100 / 0.1
B: MeOH / AcOH = 100 / 0.1
50-100%B (0-5 min), 100%B (5-6 min)
0.4 mL/min (7 MPa), 37°C
20 μL (50ng/mL
MeOH / water / AcOH = 50 / 50 / 0.1)
Single Quad. MS (ESI)