

Cadenza CD-C18


Unison UK-C8

50 x 3 mm

Application

LC-MS: Taxifolin


タキシフォリン

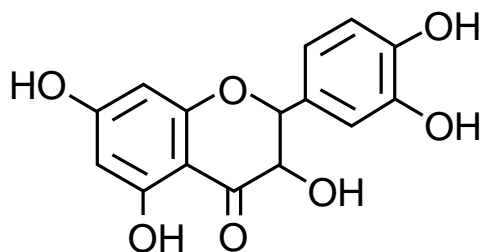

taxifolin


Cadenza CD-C18

10-100%B (0-1.5 min), 100%B (1.5-3 min)

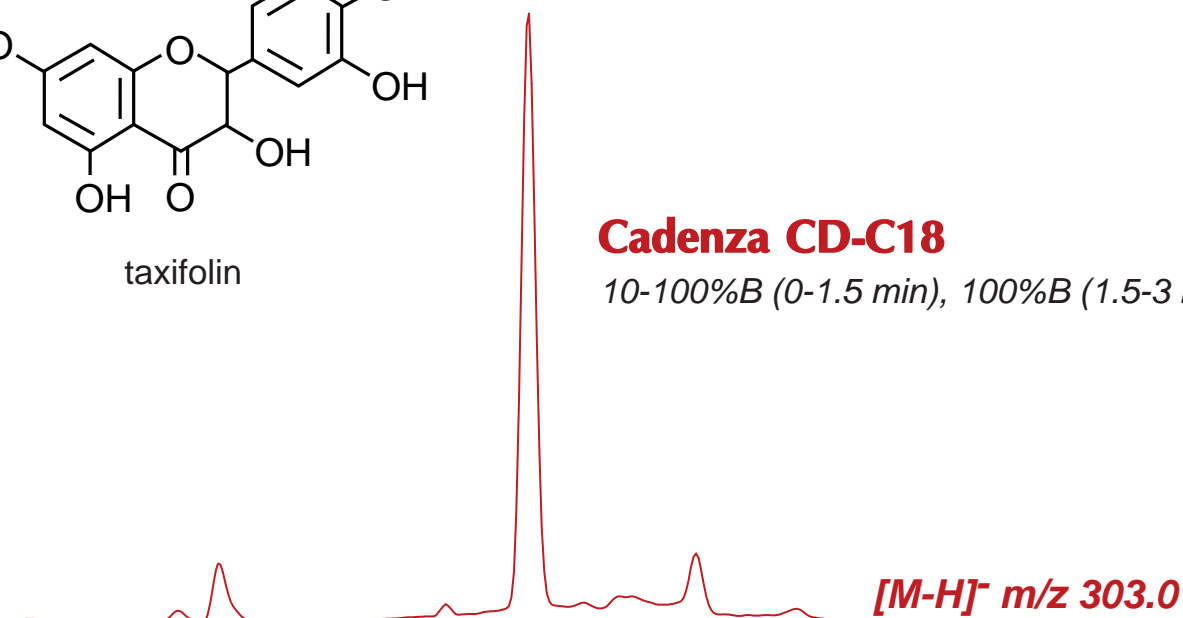

Unison UK-C8

20-100%B (0-1.5 min), 100%B (1.5-3 min)

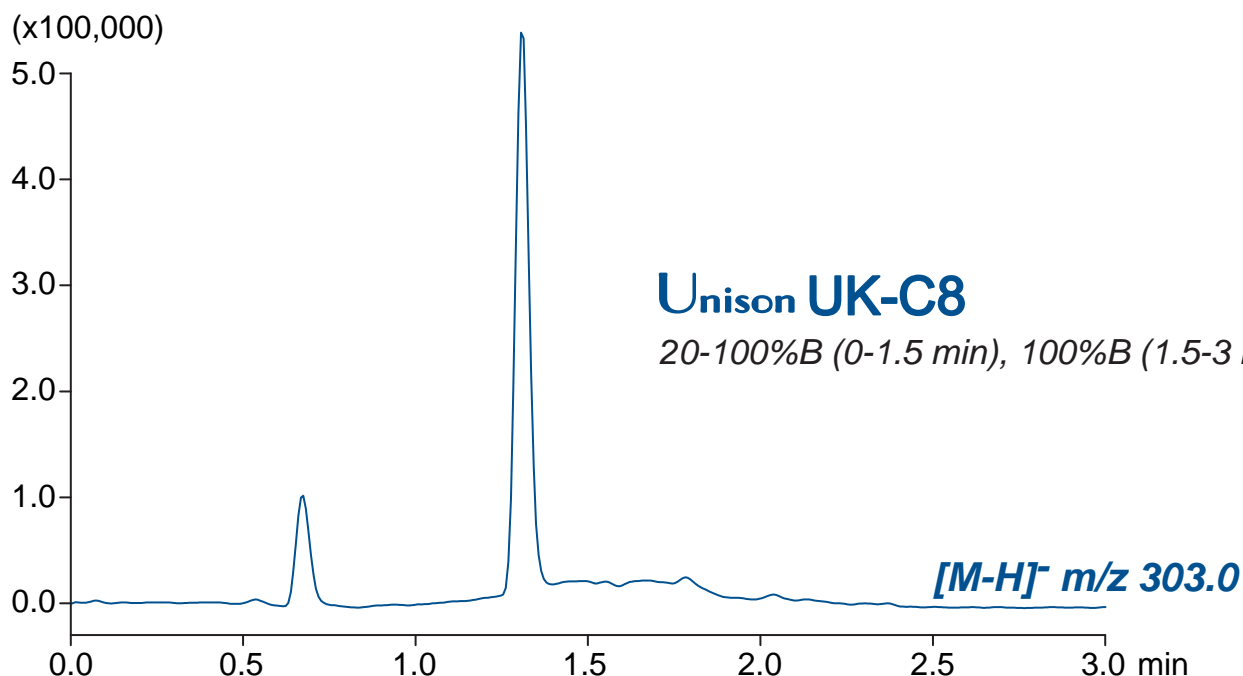

50 x 3 mm

A: water / acetic acid = 100 / 1

B: methanol / acetic acid = 100 / 1

0.4 mL/min (3-4 MPa), 50°C

10 μ L (500ng/mL, MeOH / water / AcOH = 50 / 50 / 0.1), Single Quad. MS (ESI)