

Intrada SEC

250 x 4.6 mm

Application


SEC: Polyvinylpyrrolidone (PVP)

ポリビニルピロリドン(PVP)のSEC分離




SEC: PVP

Intrada SEC

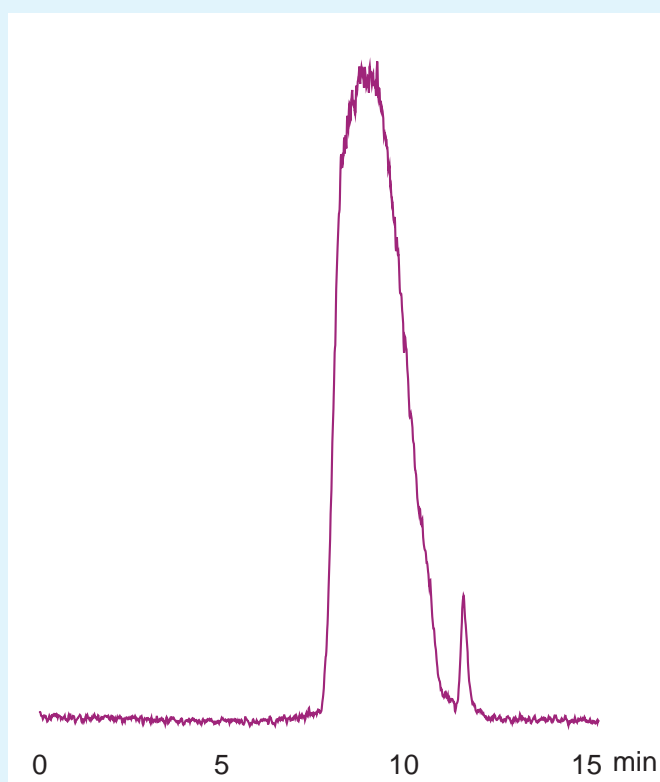

Polyvinylpyrrolidone K30

MW 40 kDa

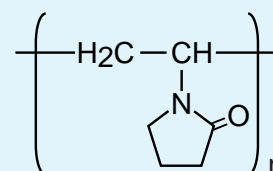

873

*Intrada SEC, 250 x 4.6 mm
 100mM HCOONH₄ /MeOH = 70 / 30
 0.3mL/min (6.3MPa), 50°C
 ELSD (spray chamber 50deg.C, drift tube 100deg.C)
 10 μL (2mg/mL, water)*