



Intrada Amino Acid

100 x 3 mm

HPLC COLUMN APPLICATION

LC-MS: Cyclic amino acids

環状アミノ酸の LC-MS 分析

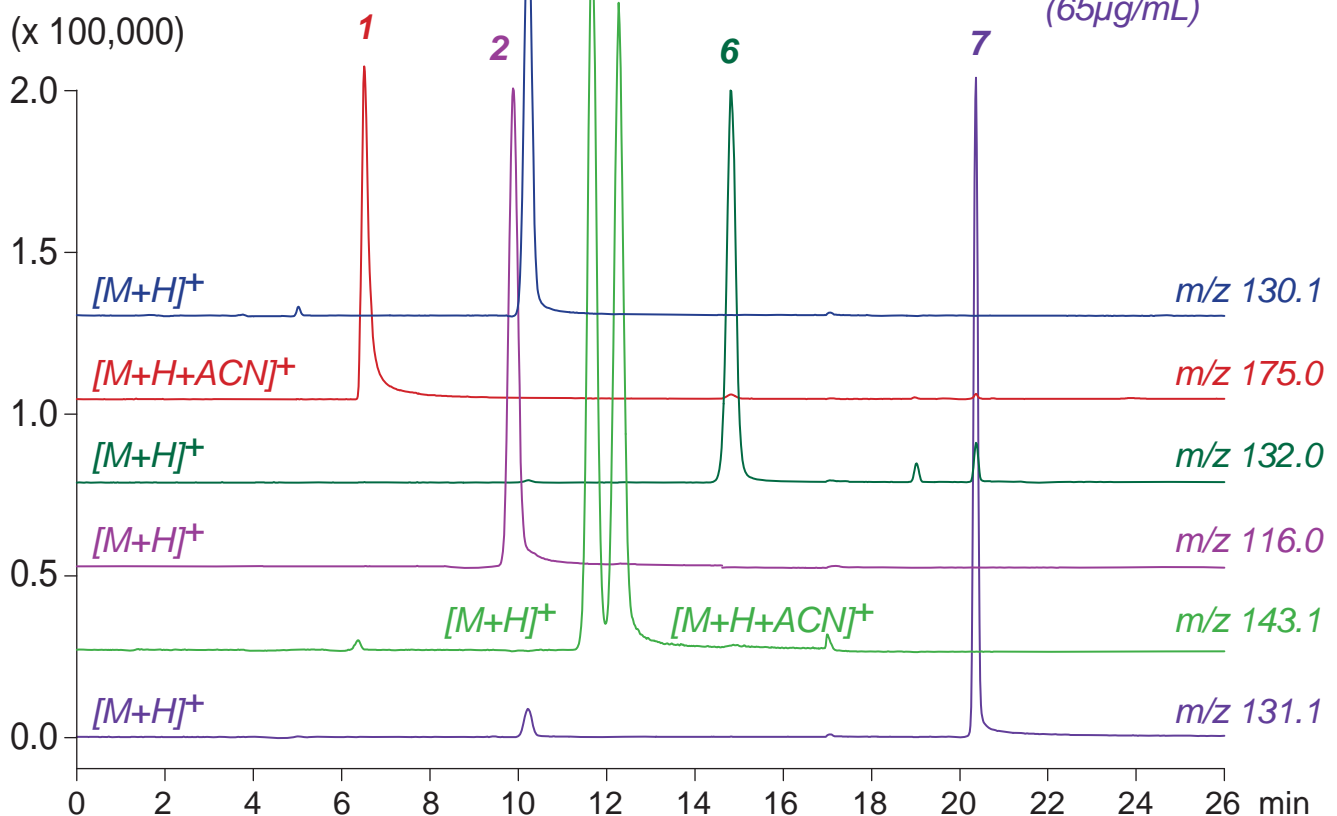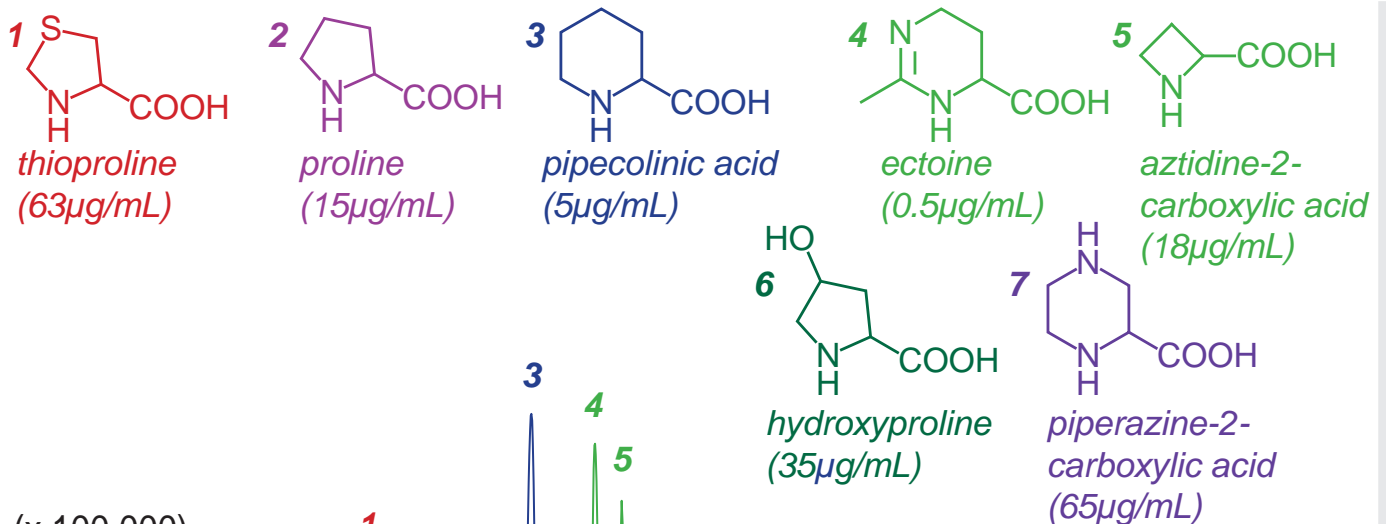

Intrada Amino Acid, 100 x 3 mm

A: acetonitrile / formic acid = 100 / 0.5

B: acetonitrile / 100mM ammonium formate = 20 / 80

15-20%B (0-15min), 95%B (15-26min)

0.4mL/min (3MPa), 37°C, Single Quad. MS (ESI)

5µL (0.5-65µg/mL, 1% HCOOH)