


Scherzo SW-C18
 Scherzo SM-C18
 Scherzo SS-C18

100 x 4.6 mm

HPLC COLUMN APPLICATION

Aminophenylboronic acid

アミノフェニルボロン酸

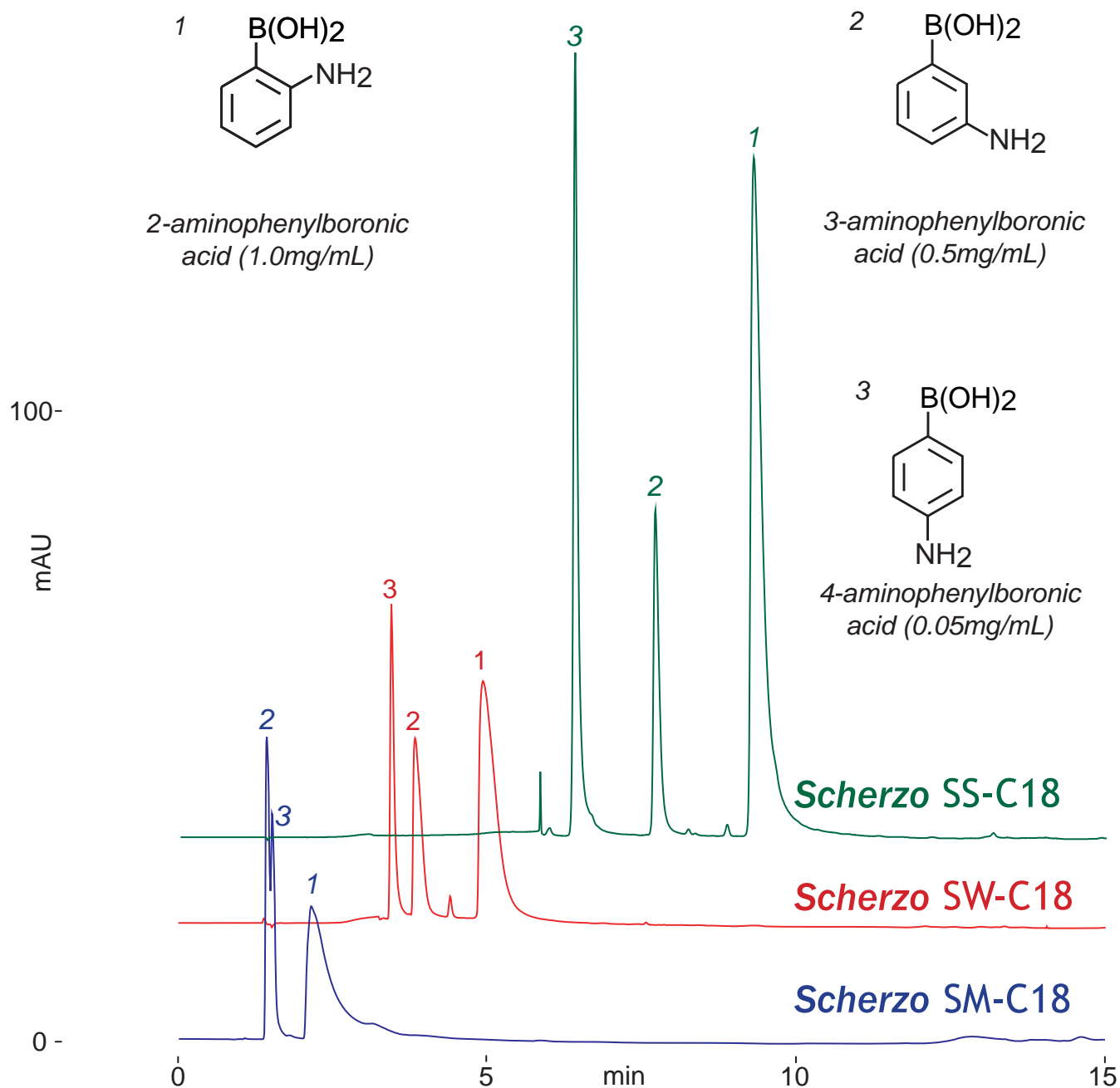

100 x 4.6 mm

A: water / formic acid = 100 / 0.1

B: 100mM ammonium formate / acetonitrile = 50 / 50

0-100%B (0-15min)

1.0mL/min (10MPa), 37°C, 260nm

10µL (0.05-0.5mg/mL, 10mM NaOH)