








Nardis ND-NX **MF**

150 x 2 mm

HPLC COLUMN APPLICATION

LC-MS: K-value (freshness index for fish) related compounds

K値 (魚鮮度指標) 関連化合物の LC-MS 分析

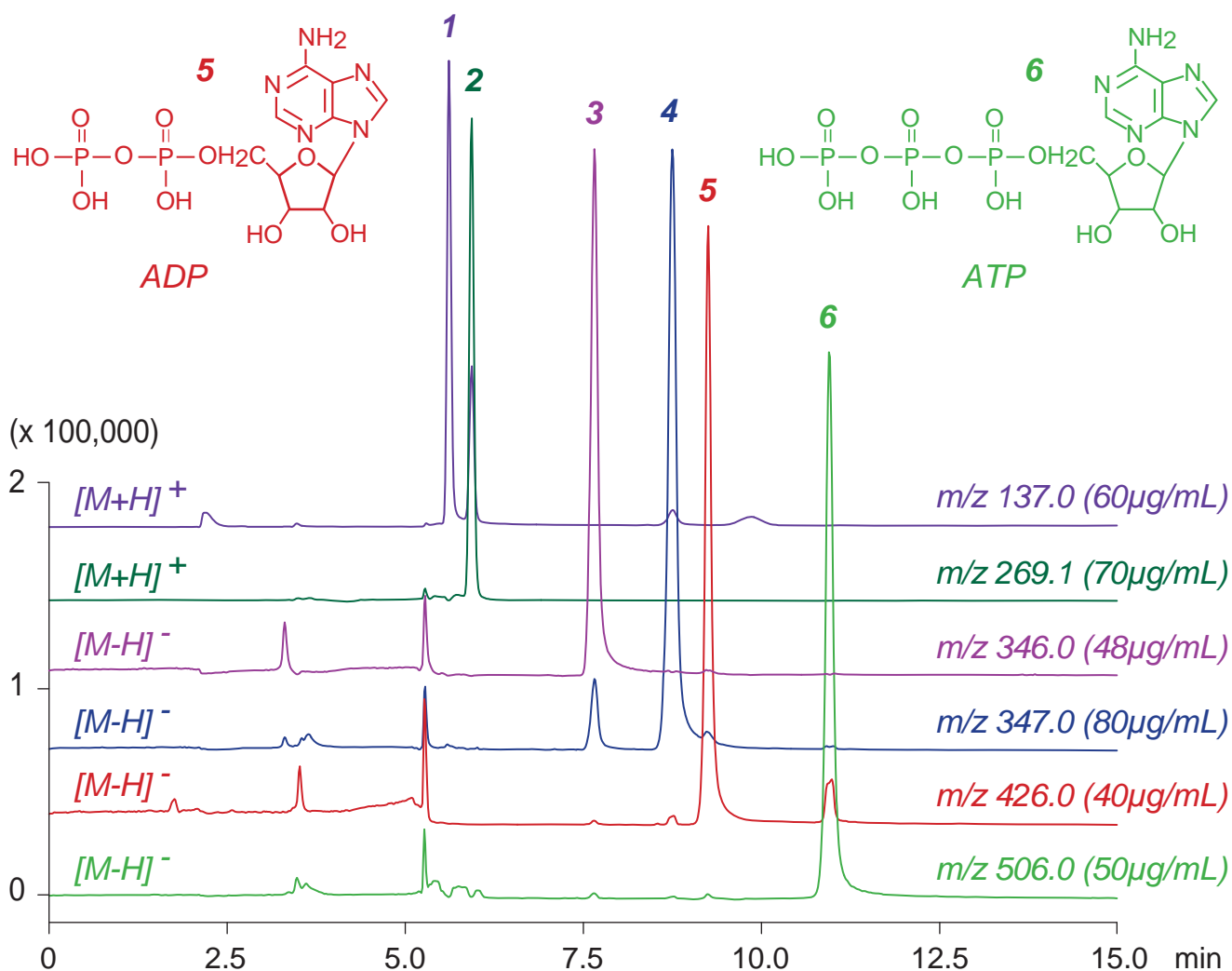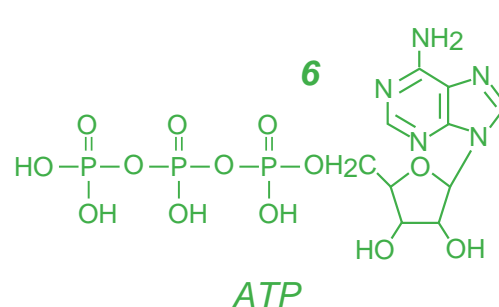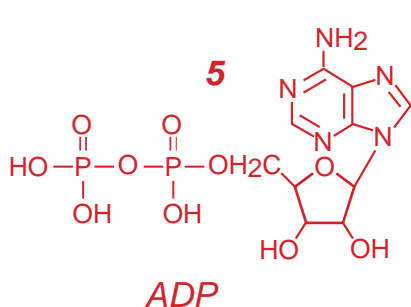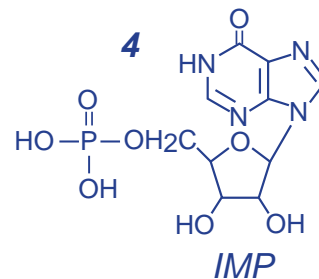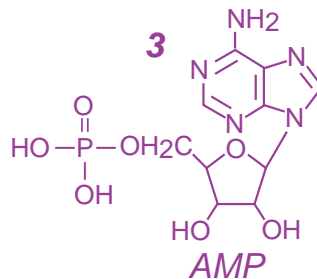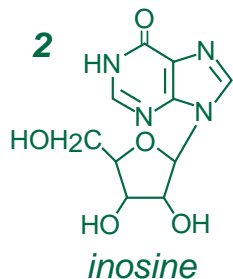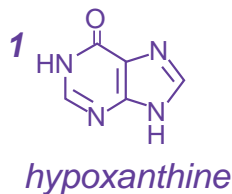

Nardis ND-NX MF, 150 x 2 mm

A: acetonitrile

B: acetonitrile / 100mM ammonium acetate = 50 / 50

5-70%B (0-2min), 70-80%B (2-15min)

0.2mL/min (4MPa), 37°C, Single Quad. MS (ESI)

2µL (40-80µg/mL, 100mM ammonium acetate)