
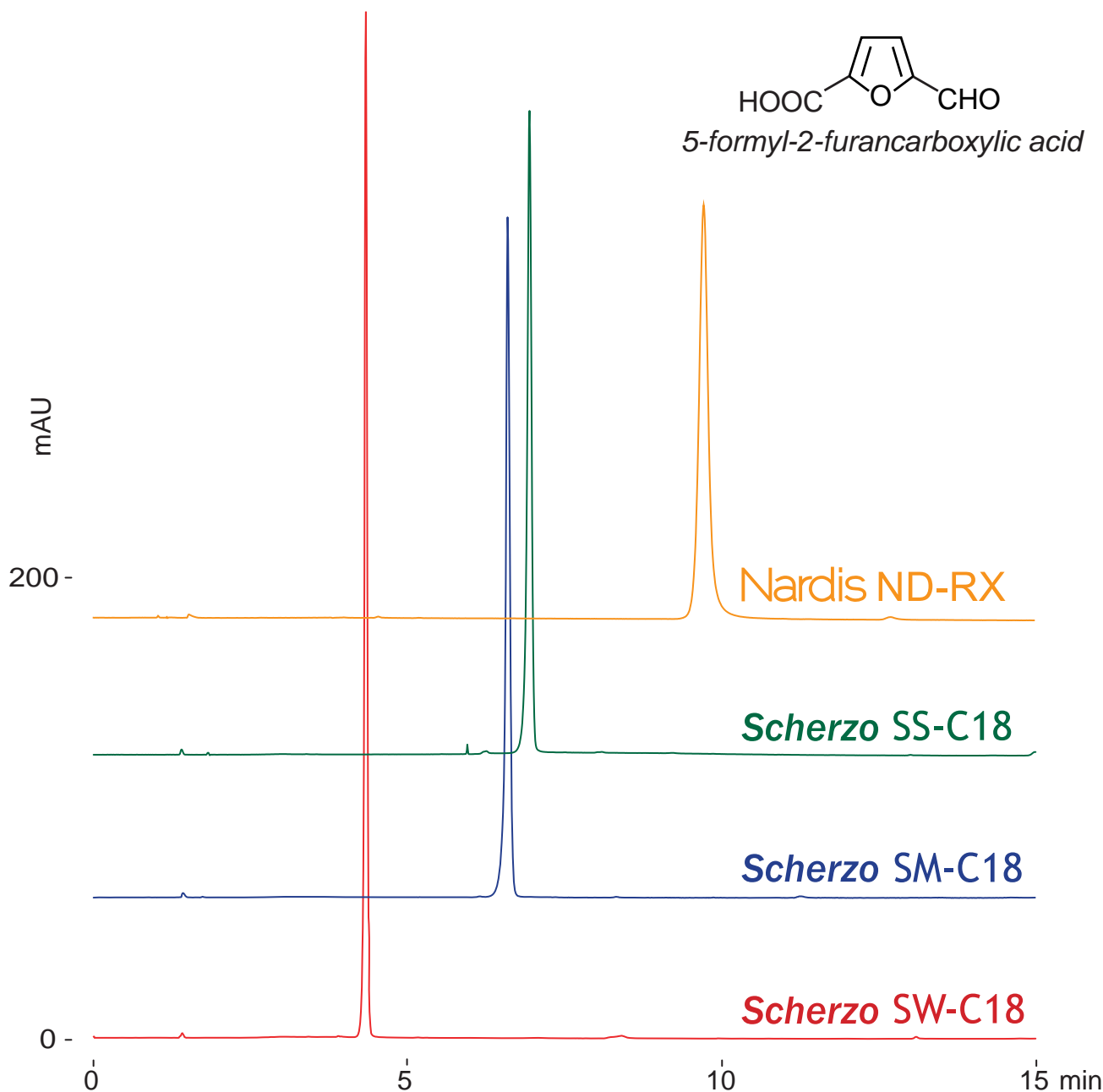


Scherzo SM-C18 Nardis ND-RX
 Scherzo SW-C18 Scherzo SS-C18 100 x 4.6 mm

HPLC COLUMN APPLICATION

5-Formyl-2-furancarboxylic acid (Furfural)
 5-ホルミル -2-フランカルボン酸 (フルフラール類)

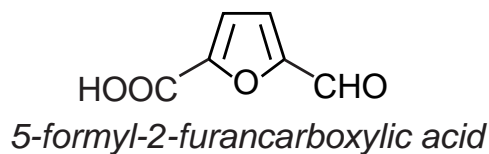

100 x 4.6 mm
 A: water / formic acid = 100 / 0.1
 B: 100mM ammonium formate / acetonitrile = 50 / 50
 0-100%B (0-15min)
 1.0 mL/min (9MPa), 37°C, 280nm
 10µL (0.035mg/mL, water / formic acid = 100 / 1)