


Cadenza CD-C18 HT

150 x 2 mm

HPLC COLUMN APPLICATION

LC-MS/MS: Antiparasitic drugs

抗寄生虫薬のLC-MS/MS分析

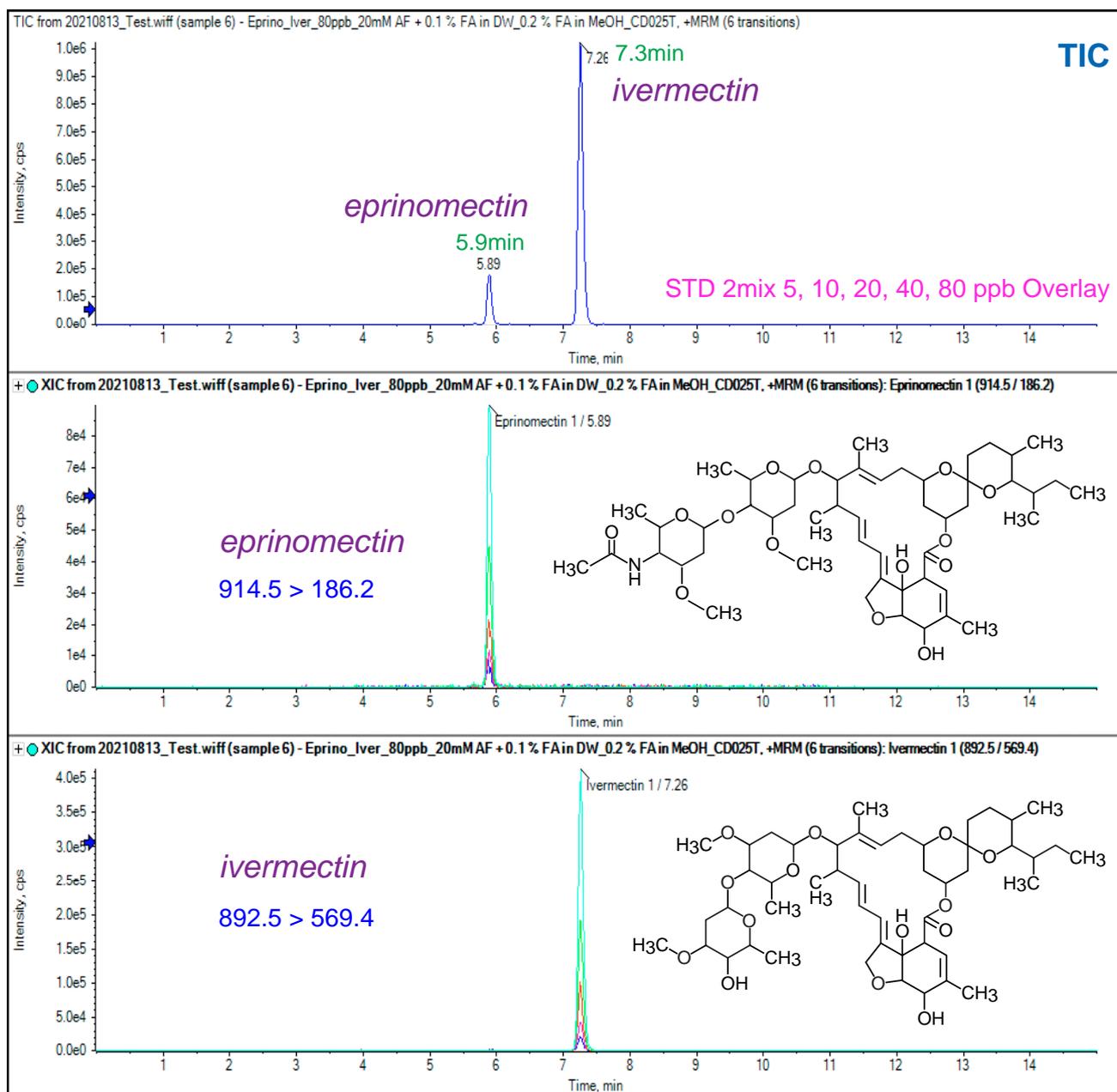

Cadenza CD-C18 HT, 150 x 2 mm

A: 0.1% formic acid + 20mM ammonium formate in water

B: 0.2% formic acid in methanol

10%B (0-0.5min), 10-95%B (0.5-2min), 95%B (2-9min),

95-10%B (9-9.1min), 10%B (9.1-15min)

0.3mL/min, 40°C, Triple Quad. MS (ESI, Positive)

5µL (5, 10, 20, 40, 80 ppb)

Courtesy of PEAKMAN SP, KOREA