


Unison UK-Phenyl Cadenza CX-C18
Unison UK-C8 Cadenza CW-C18
Unison UK-C18 Cadenza CD-C18

100 x 4.6 mm


HPLC COLUMN APPLICATION

Tris (2,2'-bipyridyl) ruthenium (II) (Dye for dye sensitized solar cell)

トリス (2,2' -ビピリジル) ルテニウム (II) (色素増感太陽電池用色素)

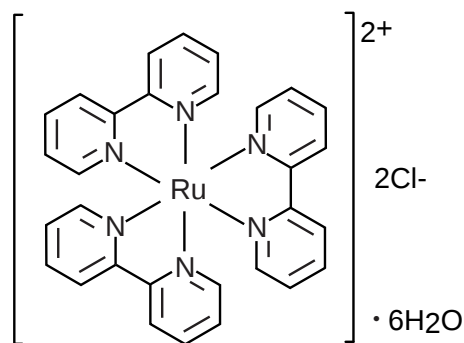

tris (2,2'-bipyridyl) ruthenium (II)
chloride hexahydrate

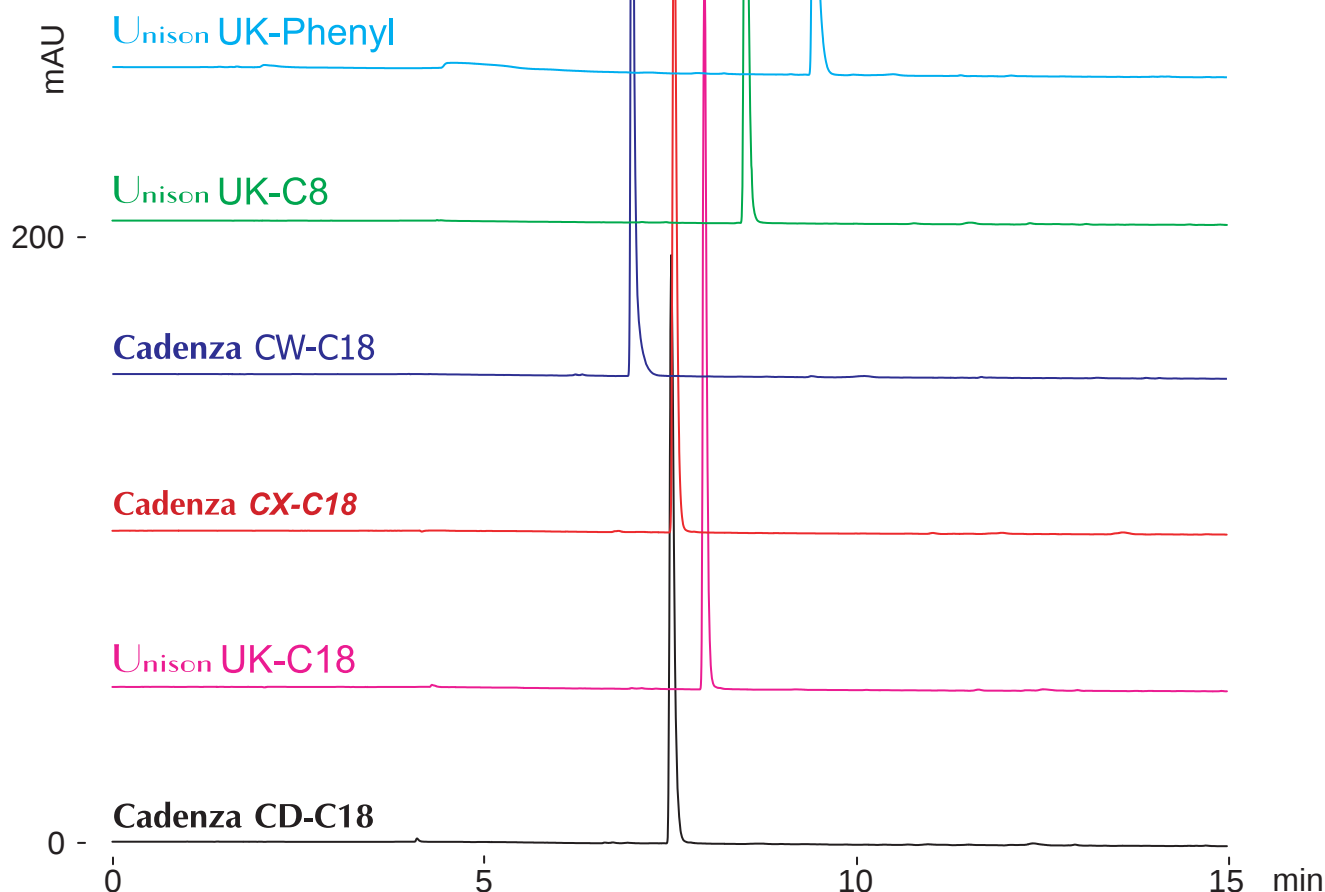

100 x 4.6 mm

A: 100mM ammonium formate

B: 100mM ammonium formate / acetonitrile = 10 / 90

0%B (0-1min), 0-60%B (1-15min)

1.0mL/min (9MPa), 37°C, 260nm

10μL (0.15mg/mL, 100mM ammonium formate)