

Scherzo SM-C18 Nardis ND-RX


Scherzo SW-C18 Scherzo SS-C18 100 x 4.6 mm

HPLC COLUMN APPLICATION


1,4-Dimethyl-1,2,4-triazolium Iodide

(Condensation organocatalysts)

1,4-ジメチル-1,2,4-トリアゾリウムヨード
(縮合有機触媒)

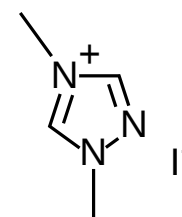

1,4-dimethyl-1,2,4-triazolium iodide

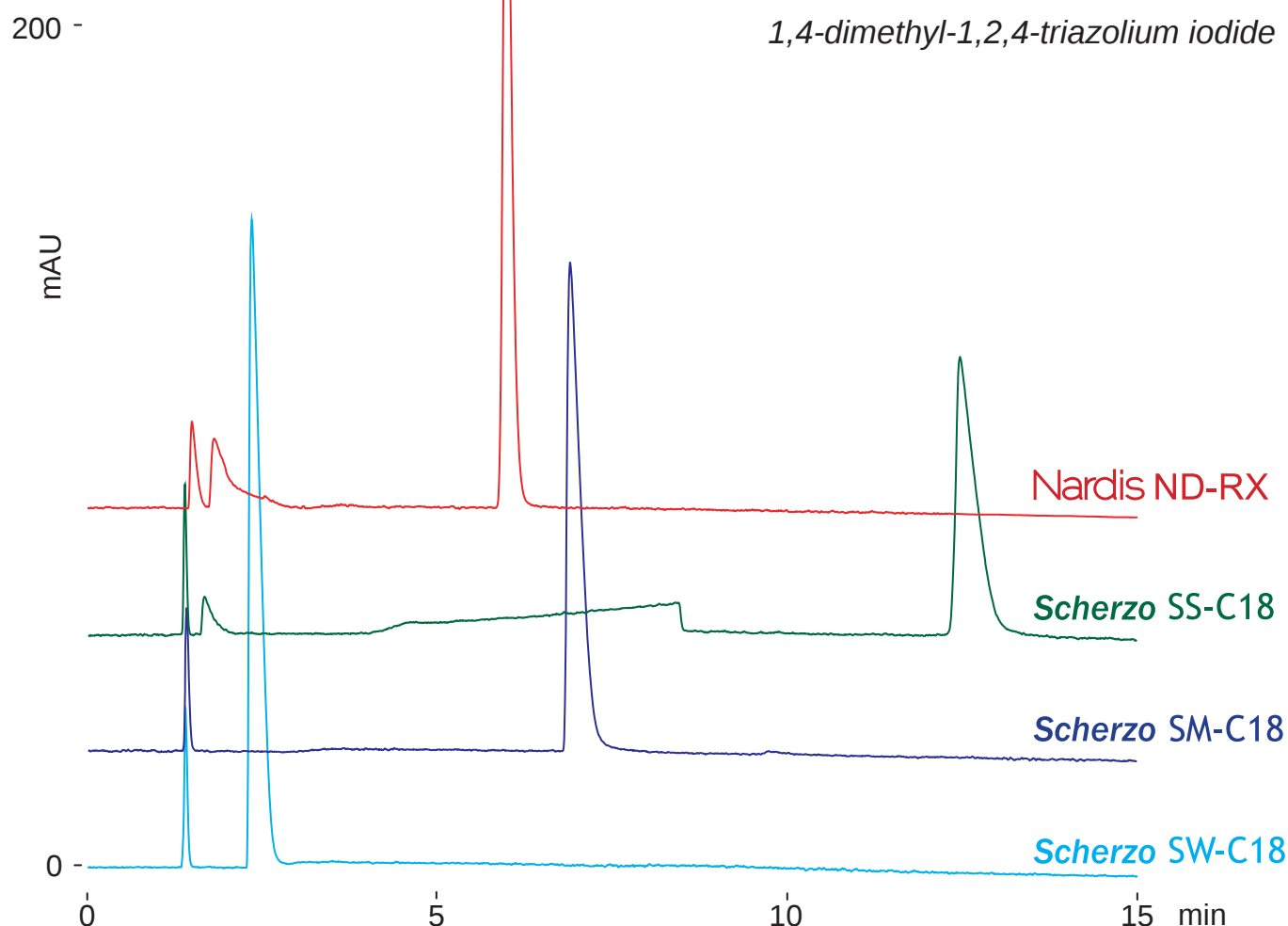

100 x 4.6 mm

A: water / formic acid = 100 / 0.1

B: 100mM ammonium formate / acetonitrile = 70 / 30

0%B (0-1min), 0-30%B (1-15min)

1.0mL/min (11MPa), 37°C, 245nm

10μL (0.5mg/mL, water / formic acid = 100 / 1)