

Cadenza CD-C18



Unison UK-C18

Unison UK-Phenyl


100 x 4.6 mm

Nitrofuran antibacterials


ニトロフラン系抗菌薬

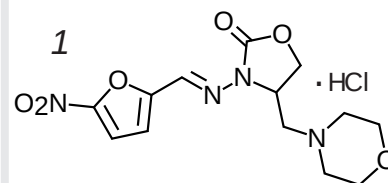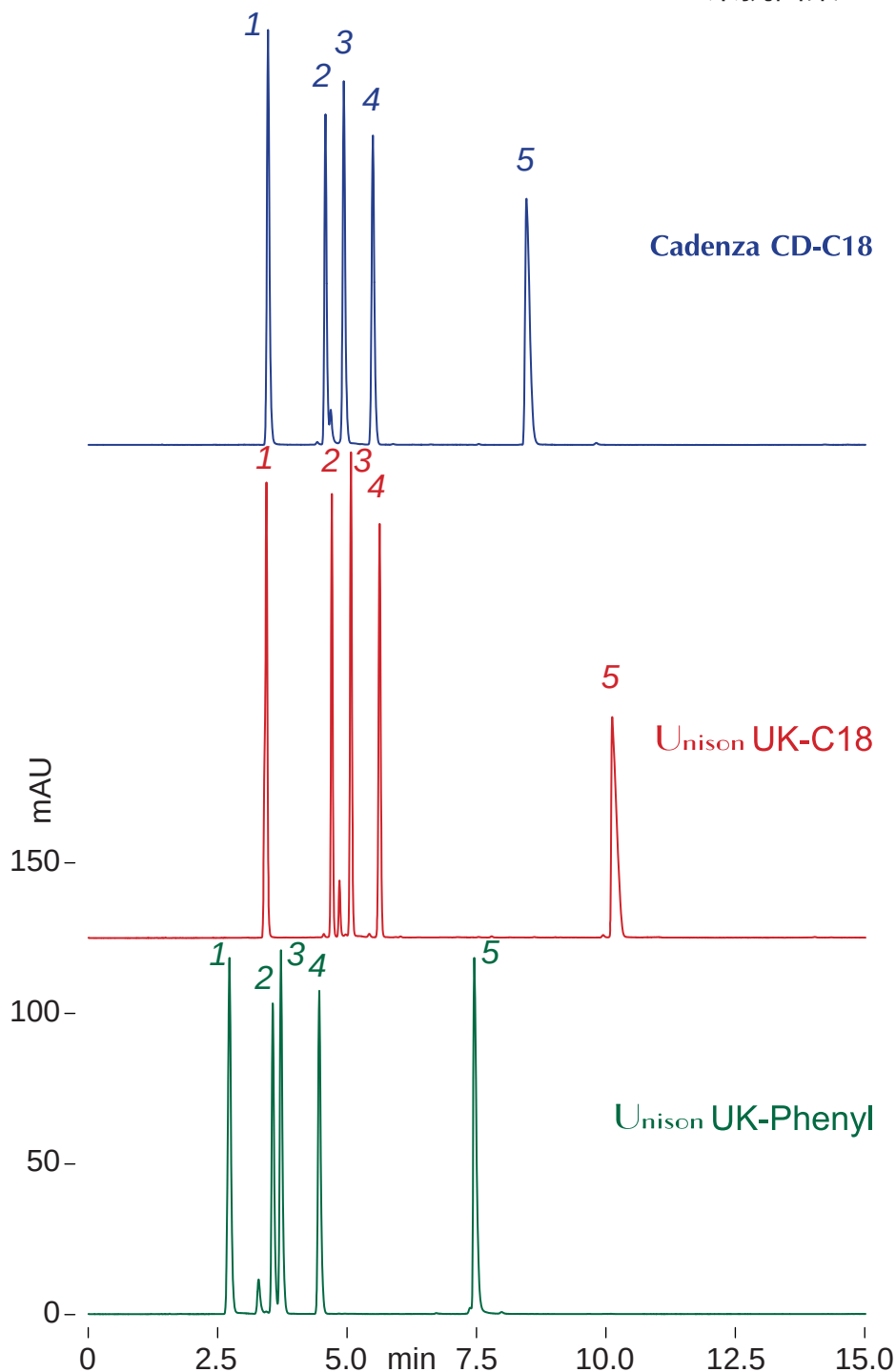


furaltadone hydrochloride
(33μg/mL)

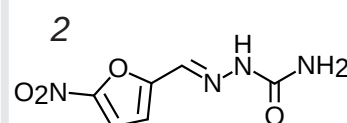


nitrofurazone (45μg/mL)

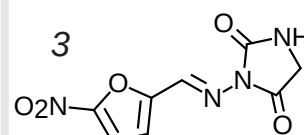

nitrofurantoin (33μg/mL)

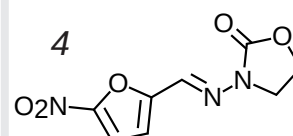

furazolidone (28μg/mL)

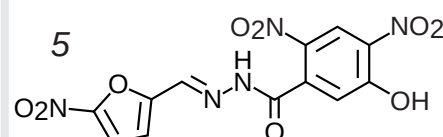

nifursol (50μg/mL)

100 x 4.6 mm

A: acetonitrile / water / formic acid = 5 / 95 / 0.1

B: acetonitrile / water / formic acid = 95 / 5 / 0.1

0-100%B (0-15min)

1mL/min (9-12MPa), 37°C, 260nm

5μL (28-50μg/mL, ACN / water / HCOOH = 40 / 60 / 1)