

Intrada Amino Acid











150 x 3 mm

Application

LC-TOFMS: Choline compounds

コリン類のLC-TOFMS分析

Metabolomics Core Facility
McGill University
GB 2017/06/16

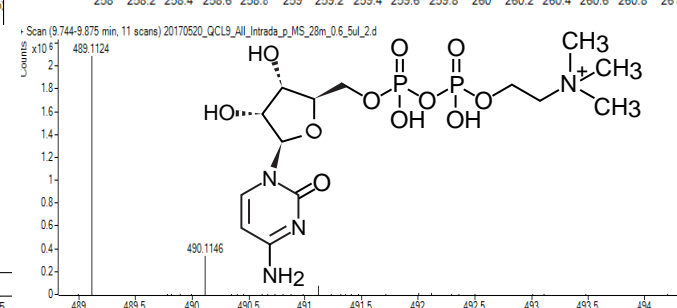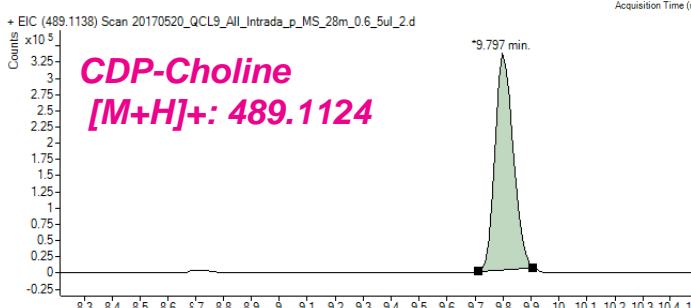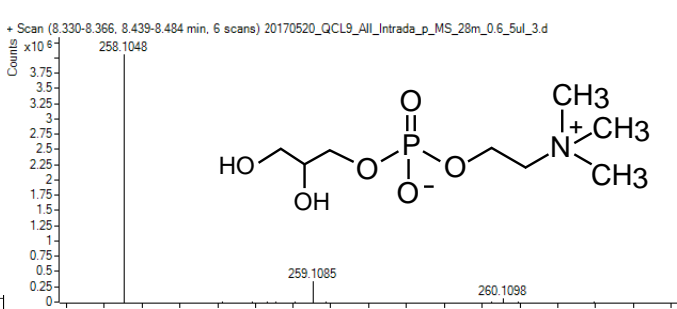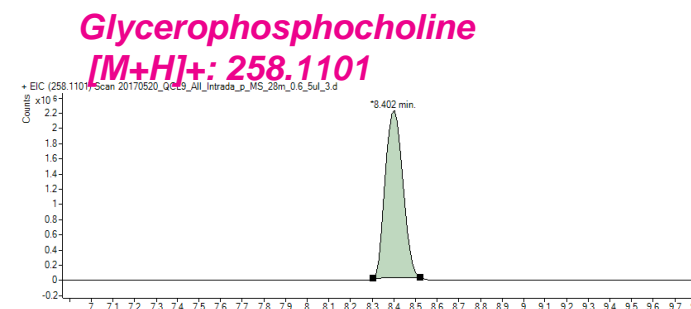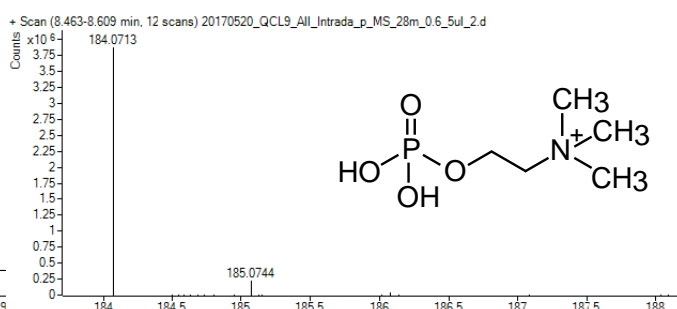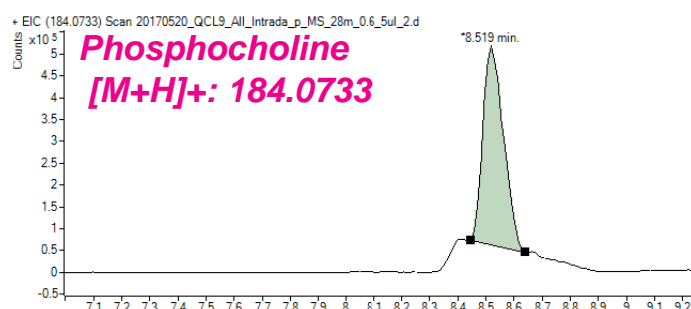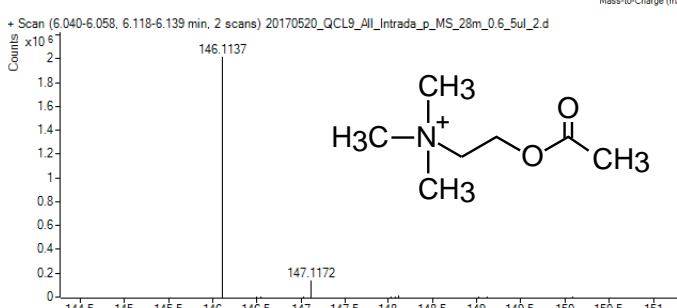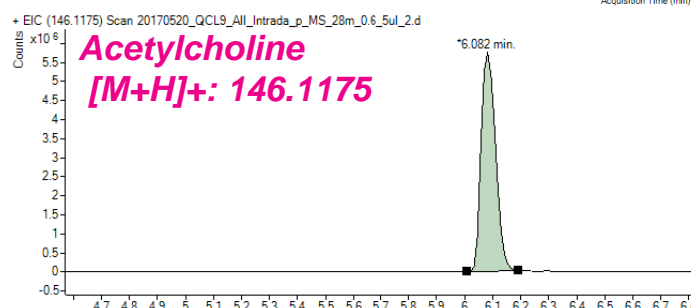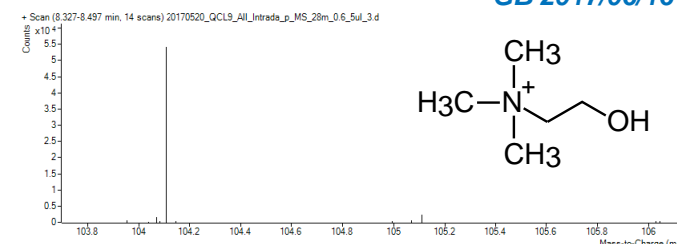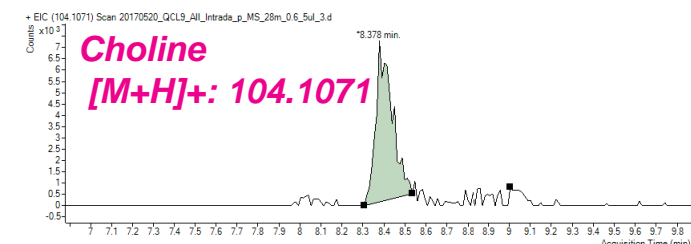

Intrada Amino Acid, 150x3mm

A) ACN / HCOOH = 100 / 0.3, B) 100mM HCOONH₄ in 20% ACN / 80% water
0-27%B (0-3min), 27-100%B (3-19.5min), 100%B (19.5-25min), 0%B (25-32min)
0.6 mL/min, 10°C, 5µL (ca. 1mM in water)

Agilent 1290 Infinity UHPLC - 6540 UHD Accurate-Mass Q-TOF mass spectrometer (ESI+)

Courtesy of Dr.Gaëlle Bridon and Dr.Daina Avizonis, McGill University, CANADA