


Nardis ND-NX

150 x 2 mm

HPLC COLUMN APPLICATION

Pyrophosphoric acid, phosphoric acid

ピロリン酸、リン酸

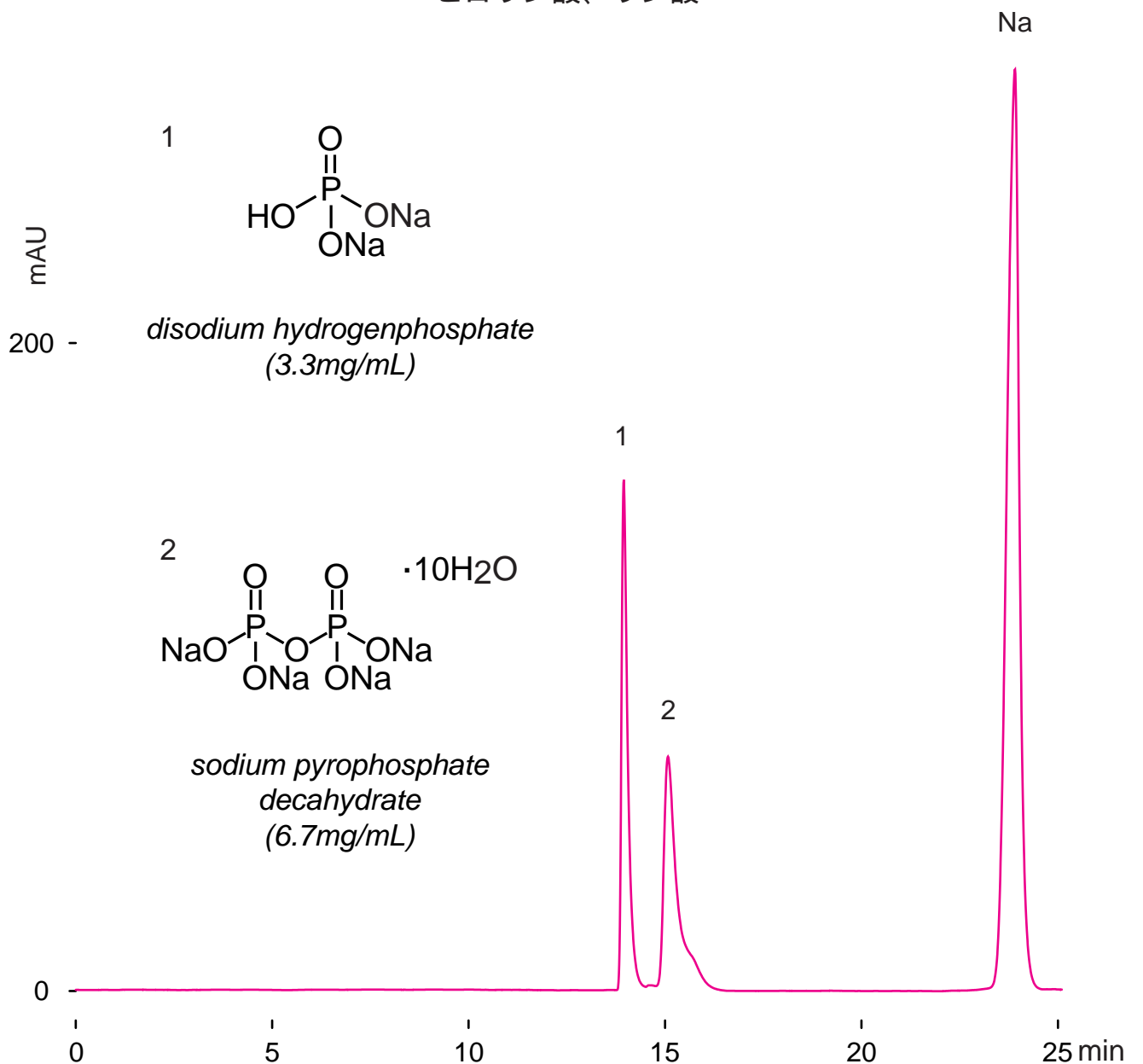

Nardis ND-NX, 150 x 2 mm

A: acetonitrile / 100mM ammonium acetate = 90 / 10

B: acetonitrile / 100mM ammonium acetate = 10 / 90

0-90%B (0-15min), 90%B (15-25min)

0.2mL/min (4MPa), 37°C,

ELSD (spray chamber 70°C, drift chamber 100°C)

3µL (3.3-6.7mg/mL, water)