

Unison UK-Phenyl Cadenza CX-C18


Unison UK-C8 Cadenza CW-C18

Unison UK-C18 Cadenza CD-C18 100 x 4.6 mm


HPLC COLUMN APPLICATION

Oxacillin (beta-Lactam antibiotic)

オキサシリン (β-ラクタム系抗生物質)

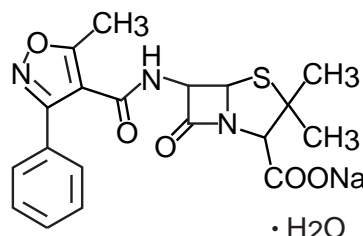

oxacillin sodium monohydrate

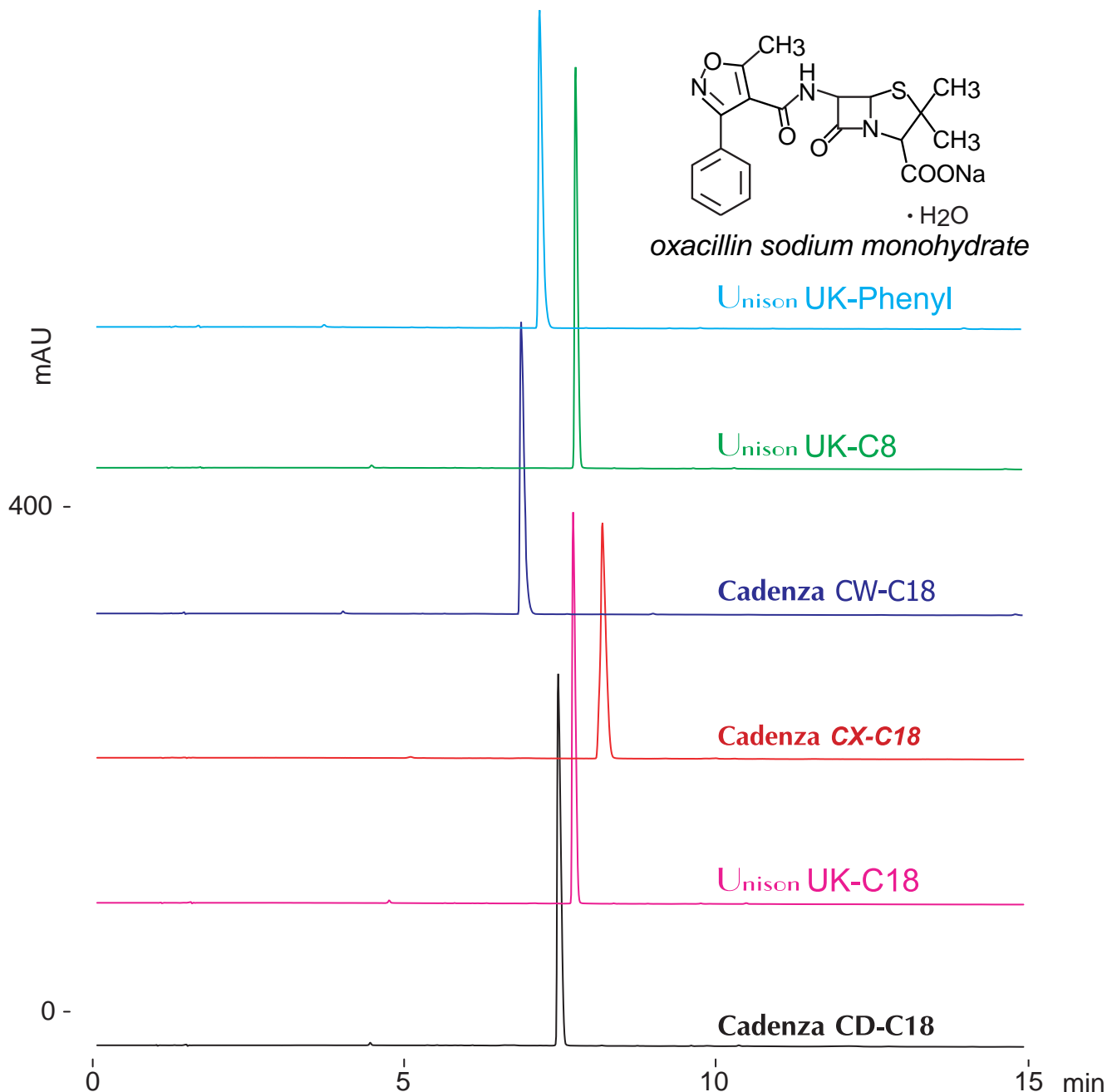

100 x 4.6 mm

A: 100mM ammonium formate / acetonitrile = 90 / 10

B: 100mM ammonium formate / acetonitrile = 10 / 90

0-70%B (0-15min)

1.0mL/min (10MPa), 37°C, 260nm

10μL (1mg/mL, 100mM ammonium formate)