

Unison UK-Phenyl Cadenza CX-C18



Unison UK-C8 Cadenza CW-C18

Unison UK-C18 Cadenza CD-C18 100 x 4.6 mm

HPLC COLUMN APPLICATION

Bergenin (Isocoumarin derivatives)

ベルゲニン (イソクマリン誘導体)

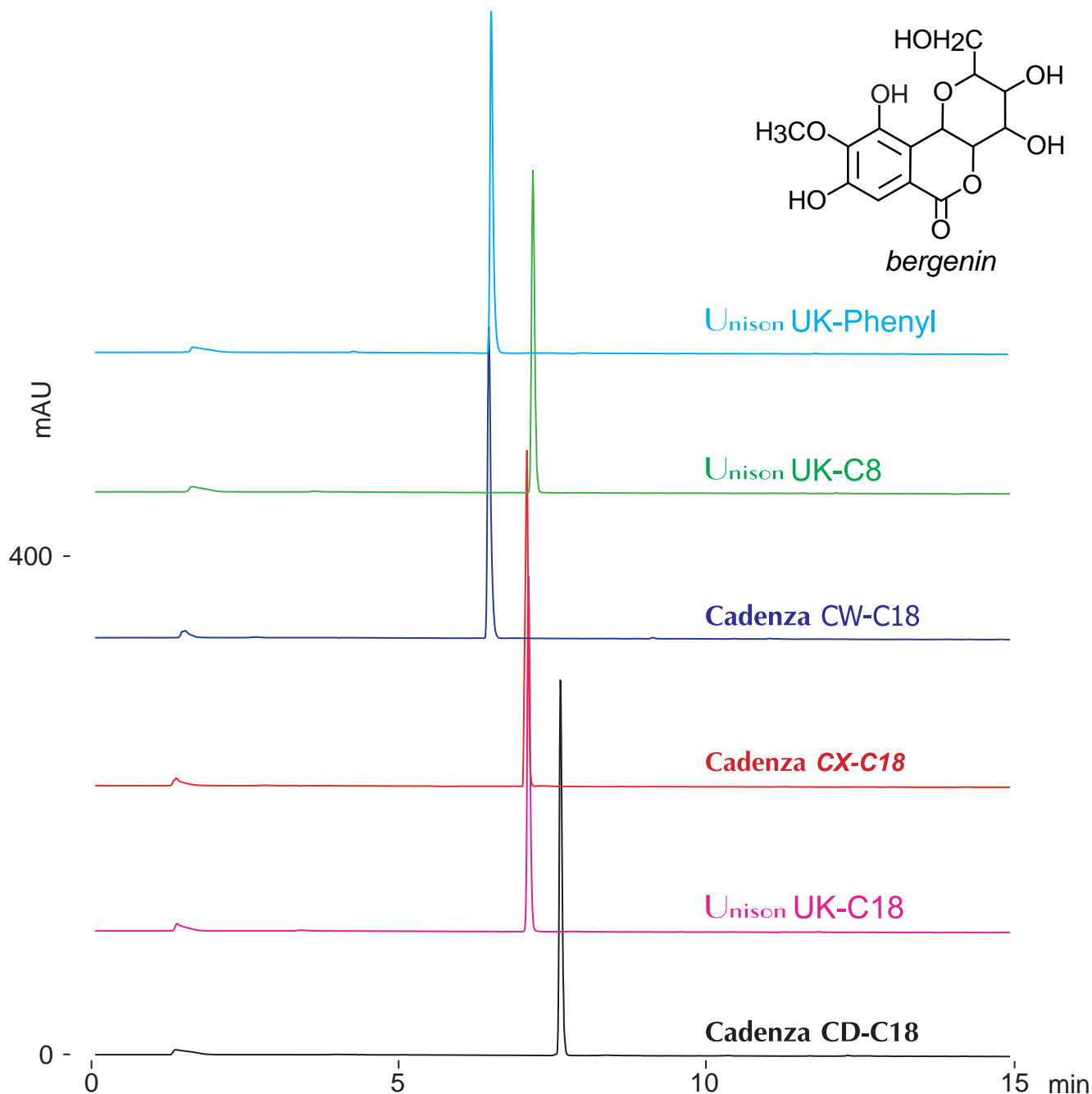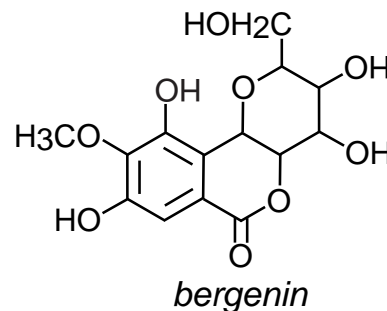

100 x 4.6 mm

A: 100mM ammonium formate

B: 100mM ammonium formate / acetonitrile = 10 / 90

0%B (0-1min), 0-40%B (1-15min)

1.0mL/min (10MPa), 37°C, 270nm

10µL (0.15mg/mL, 100mM ammonium formate / dimethyl sulfoxide = 70 / 30)