






Nardis ND-NX

150 x 3 mm

HPLC COLUMN APPLICATION

LC-MS: Aminoglycoside antibiotics

アミノグリコシド系抗生物質の LC-MS 分析

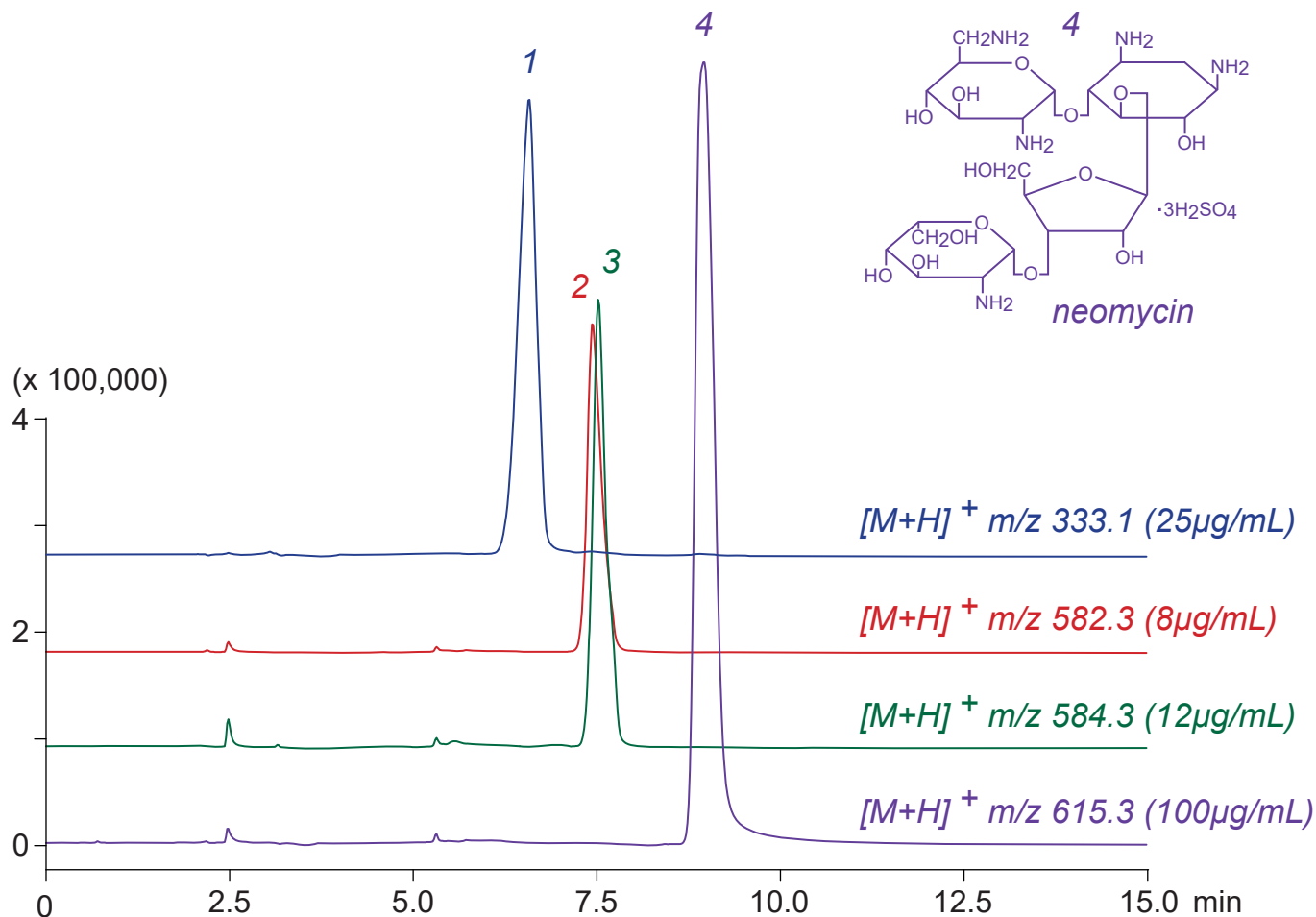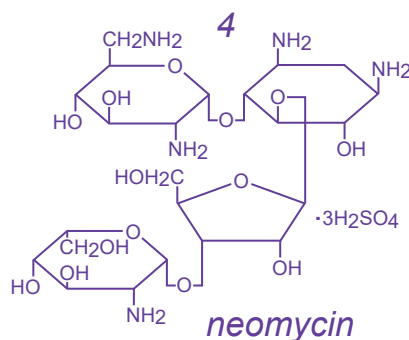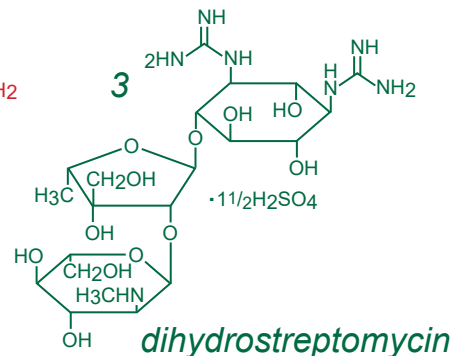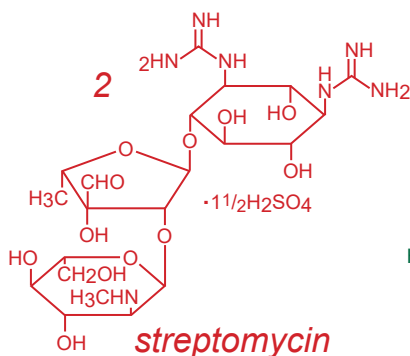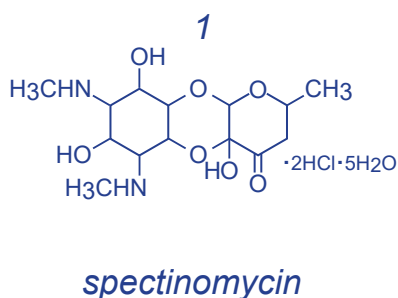

Nardis ND-NX, 150 x 3 mm

A: acetonitrile / formic acid = 100 / 0.5

B: acetonitrile / water / formic acid = 50 / 50 / 0.5

0%B (0-1min), 0-100%B (1-10min), 100%B (10-15min)

0.4mL/min (4MPa), 37°C, Single Quad. MS (ESI)

10 μ L (8-100 μ g/mL, 0.1% formic acid)