

**Intrada Organic Acid**



**Nardis ND-NX MF**

150 x 2 mm


**HPLC COLUMN APPLICATION**

**LC-MS: Coenzyme A and acetyl coenzyme A**


コエンザイム A 及びアセチルコエンザイム A (補酵素) の LC-MS 分析

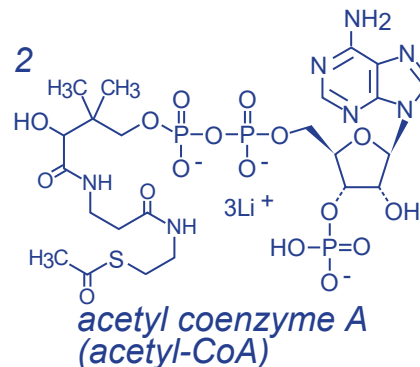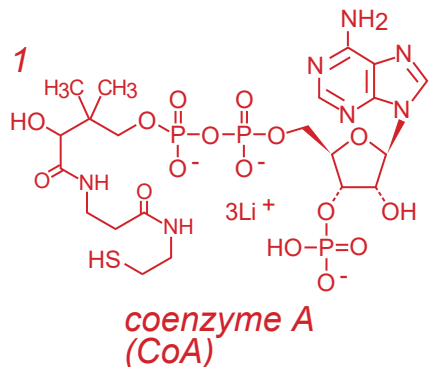

**Intrada Organic Acid**

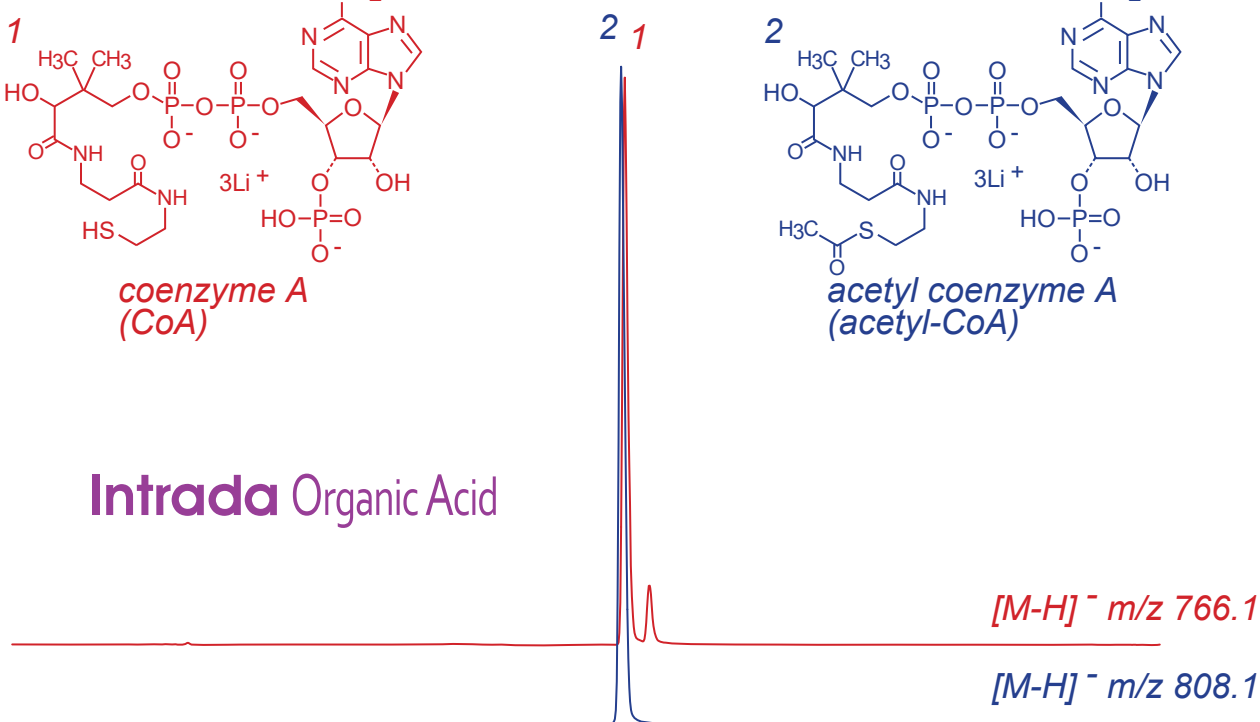

**Nardis ND-NX MF**

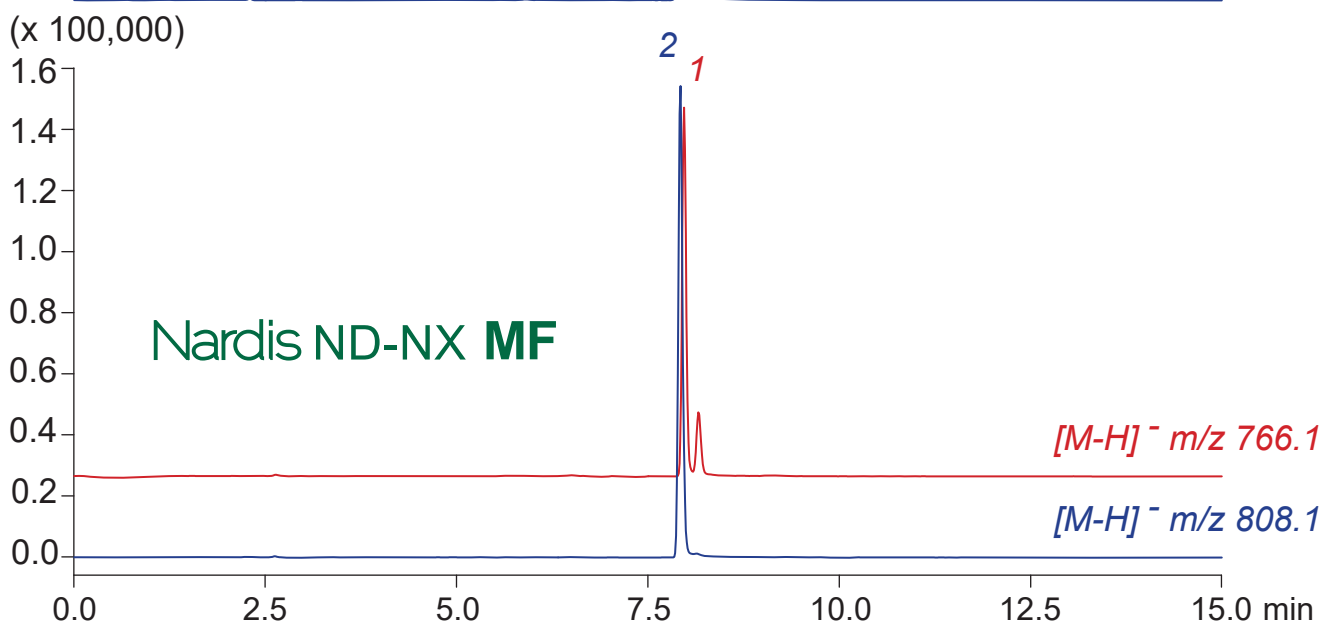

150 x 2 mm

A: ACN / 100mM HCOONH<sub>4</sub> = 95 / 5

B: ACN / 100mM HCOONH<sub>4</sub> = 10 / 90

0%B (0-1min), 0-80%B (1-7min), 80%B (7-15min)

0.2mL/min (4-5MPa), 37°C, Single Quad. MS (ESI)

5μL (20μg/mL, 50mM ammonium formate)