


Unison UK-C18 HT

150 x 2 mm

HPLC COLUMN APPLICATION

LC-MS/MS: Chlorophenols

クロロフェノール類の LC-MS/MS 分析

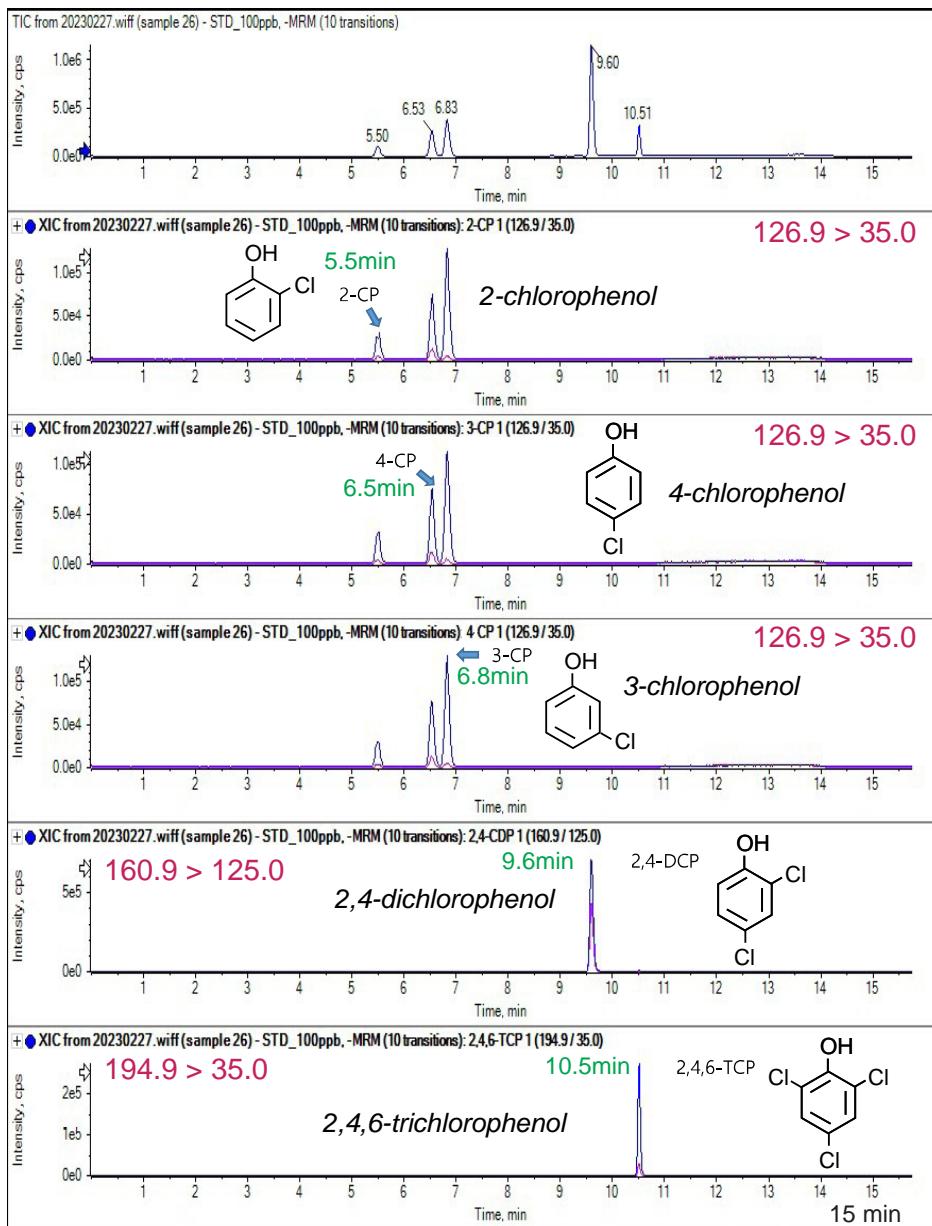

Unison UK-C18 HT, 150 x 2 mm

A: 5mM ammonium formate

B: methanol

40-55% B (0-6min), 55-70% B (6-8min), 70-95% B (8-9min),
 95% B (9-12min), 95-40% B (12-12.1min), 40% B (12.1-17min)
 0.2mL/min, 40°C, Triple Quad. MS (APCI, Negative)
 5µL (100ppb)

Courtesy of PEAKMAN SP, KOREA