

Scherzo SM-C18 MF


100 x 2 mm

HPLC COLUMN APPLICATION

LC-MS/MS: Glyphosate

グリホサートの LC-MS/MS 分析

Glyphosate STD 10 ppb

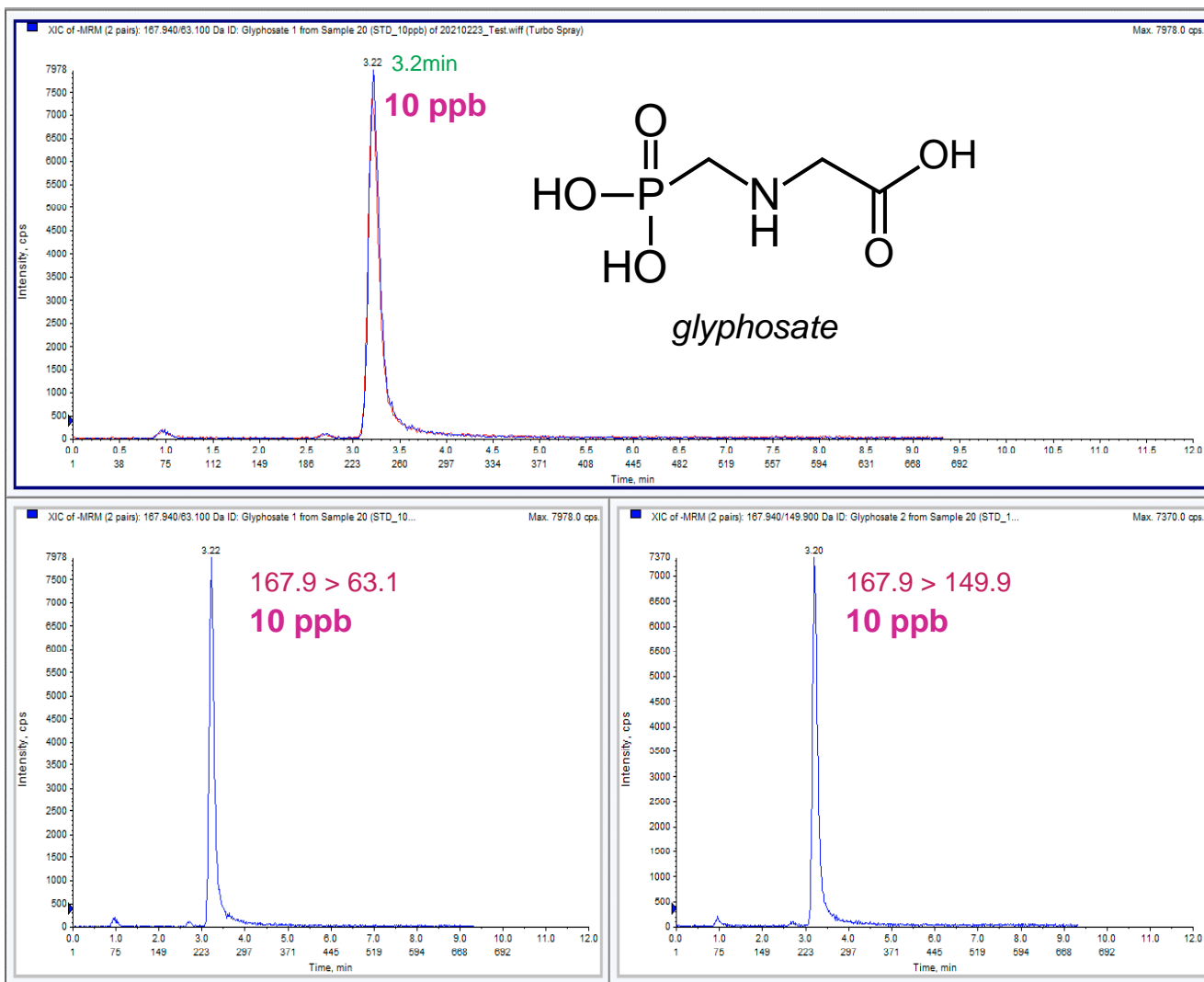

Scherzo SM-C18 MF, 100 x 2 mm

A: 0.1% formic acid in 20% methanol

B: 1% formic acid in 20% methanol

0%B (0-1min), 0-100%B (1-4min), 100%B (4-8min),

100-0%B (8-8.1min), 0%B (8.1-15min)

0.25mL/min, 40°C, Triple Quad. MS (ESI, Negative)

5μL (10 ppb)

Courtesy of PEAKMAN SP, KOREA