


Cadenza CD-C18 HT

150 x 2 mm

HPLC COLUMN APPLICATION

LC-MS/MS: Aflatoxin
アフラトキシンの LC-MS/MS 分析

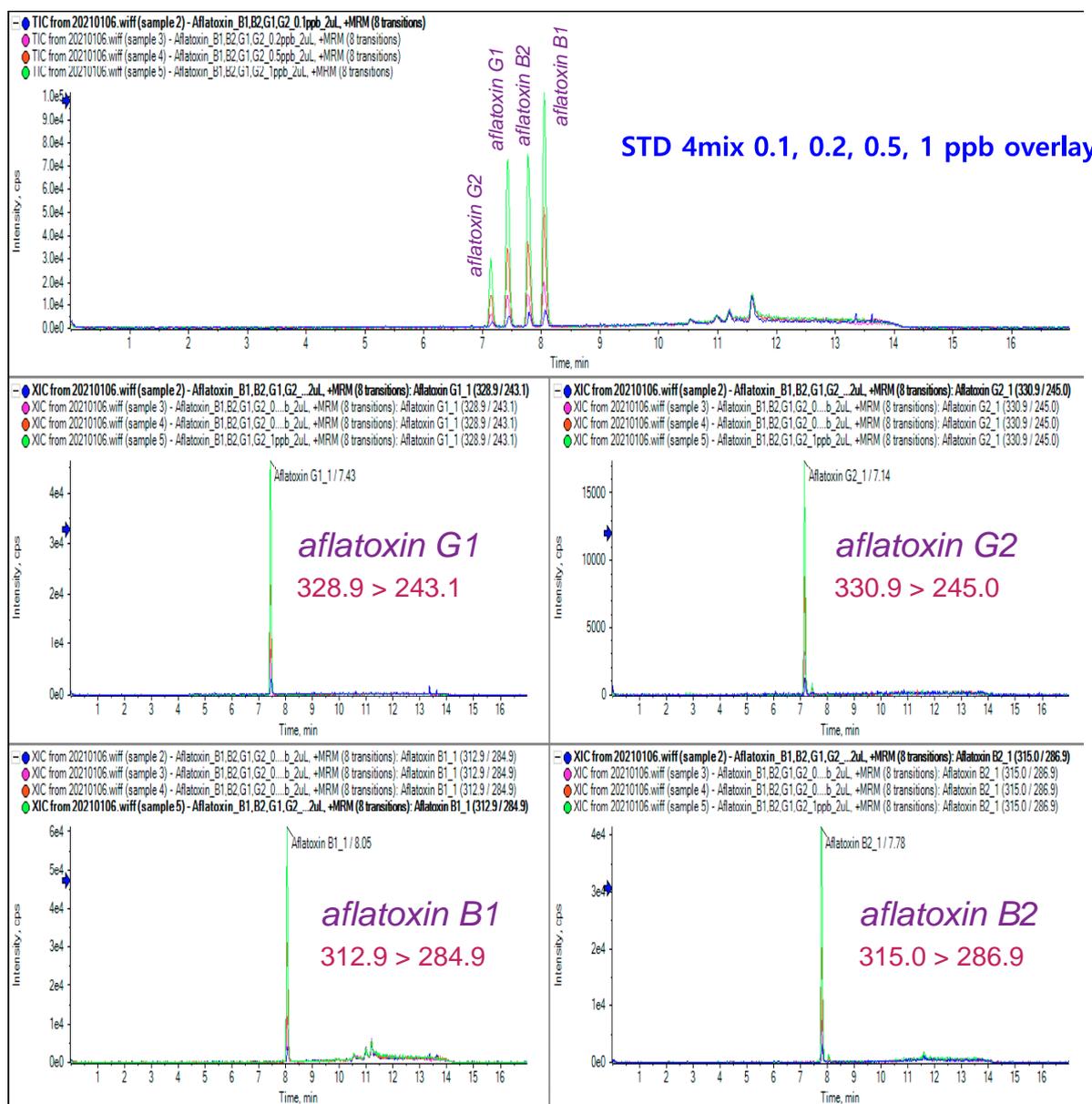

Cadenza CD-C18, 150 x 2 mm

A: 0.1% formic acid + 5mM ammonium acetate in water

B: 0.1% formic acid + 5mM ammonium acetate in methanol

5%B (0-0.5min), 5-40%B (0.5-2min), 40-100%B (2-9min),

100%B (9-12min), 100-5%B (12-12.1min), 5%B (12.1-18min)

0.2mL/min, 40°C, Triple Quad. MS (ESI, Positive)

2µL (0.1, 0.2, 0.5, 1 ppb)

Courtesy of PEAKMAN SP, KOREA