


Unison UK-C18 HT

100 x 2 mm

HPLC COLUMN APPLICATION

LC-MS/MS: Mancozeb

マンゼブの LC-MS/MS 分析

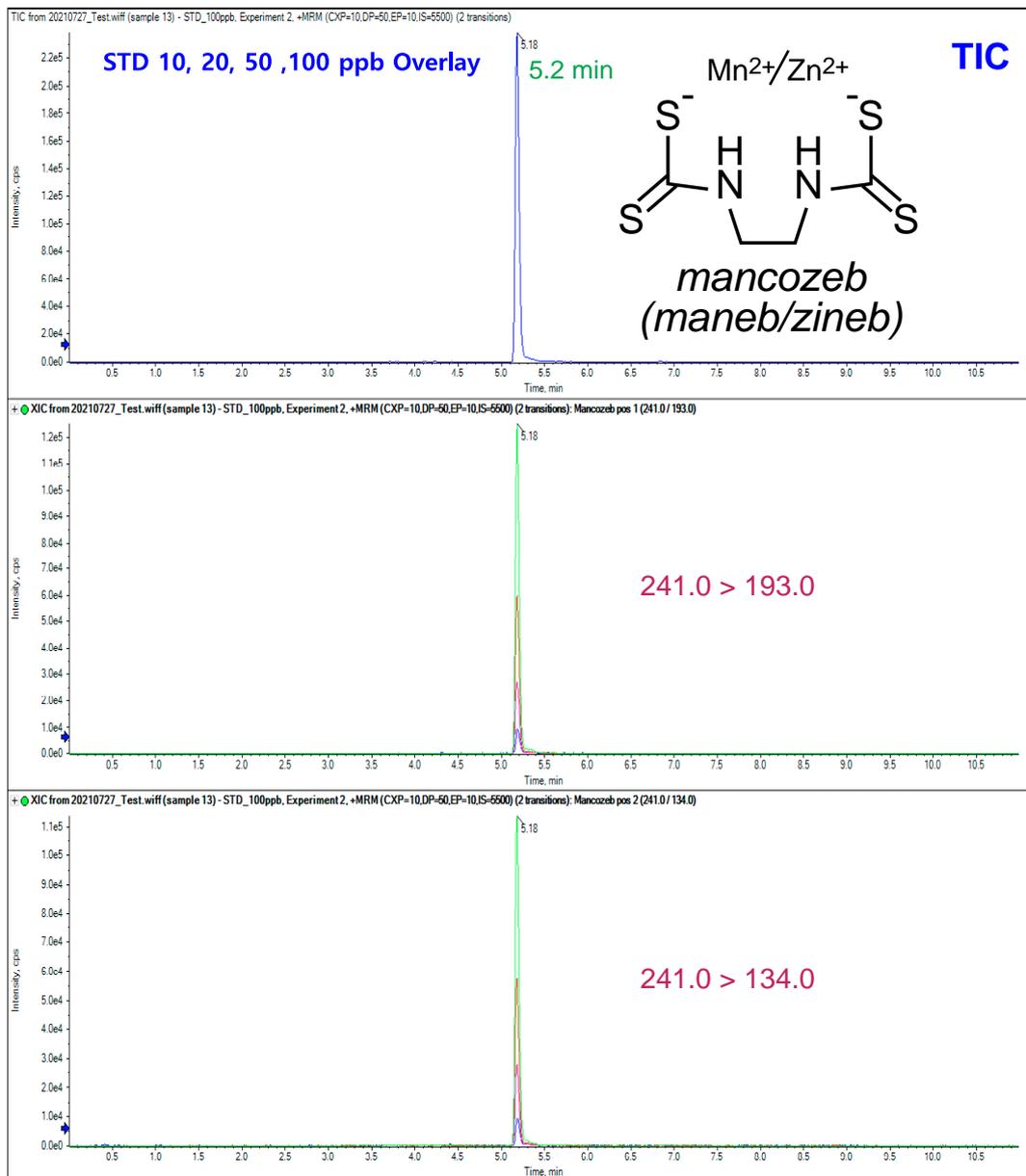

Unison UK-C18 HT, 100 x 2 mm

A: 0.1% formic acid + 5mM ammonium formate in water

B: 0.1% formic acid + 5mM ammonium formate in methanol

10%B (0-1min), 10-90%B (1-4min), 90%B (4-7min),

90-10%B (7-7.1min), 10%B (7.1-11min)

0.3mL/min, 40°C, Triple Quad. MS (ESI, Positive)

5µL (10, 20, 50, 100 ppb)

Courtesy of PEAKMAN SP, KOREA