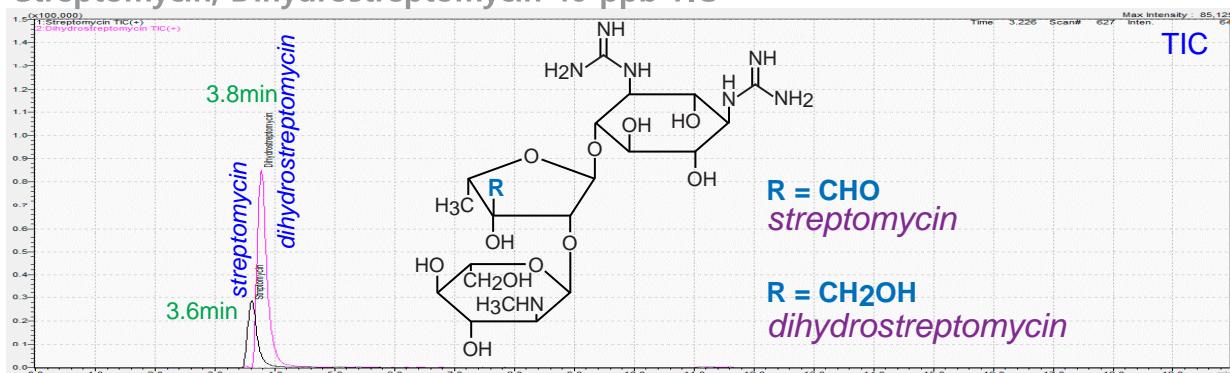
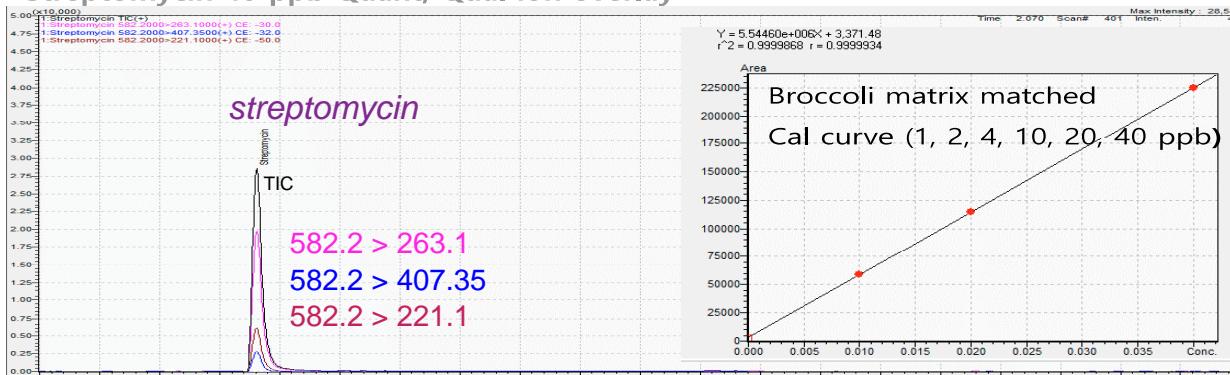
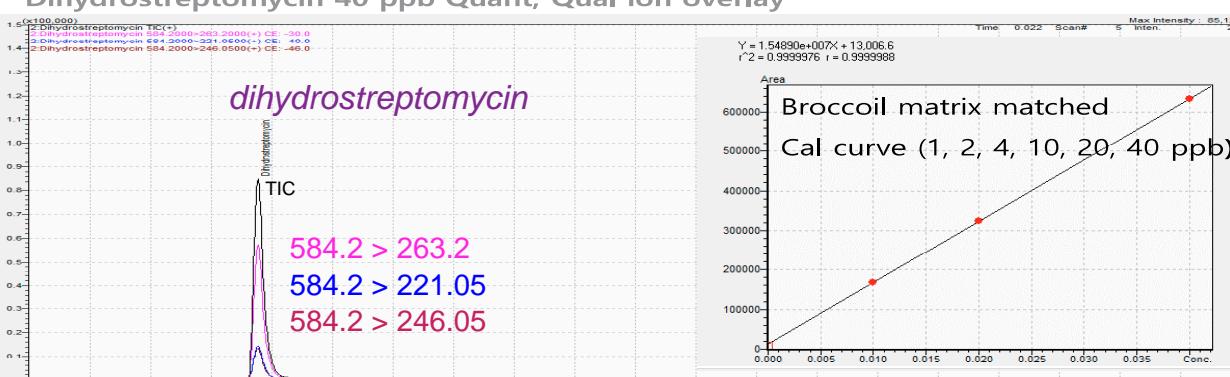


Scherzo SM-C18 MF

100 x 2 mm

HPLC COLUMN APPLICATION**LC-MS/MS: Streptomycin, dihydrostreptomycin**

ストレプトマイシン、ジヒドロストレプトマイシンの LC-MS/MS 分析

Streptomycin, Dihydrostreptomycin 40 ppb TIC**Streptomycin 40 ppb Quant, Qual ion overlay****Dihydrostreptomycin 40 ppb Quant, Qual ion overlay****Scherzo SM-C18 MF, 100 x 2 mm**

A: 30mM ammonium acetate

B: 0.8% acetic acid in 20% acetonitrile

5-70% B (0-8min), 70-100% B (8-10min), 100% B (10-13min),

100-5% B (13-13.1min), 5% B (13.1-20min)

0.2mL/min, 37°C, Triple Quad. MS (ESI, Positive)

2 μ L (40 ppb)

Courtesy of PEAKMAN SP, KOREA